

CONEX-PSD

**Two-Axis Position
& Power Sensing device**



 **Newport®**

**Command Interface
Manual**

V3.0.x

©2017 by Newport Corporation, Irvine, CA. All rights reserved.

Original instructions.

No part of this document may be reproduced or copied without the prior written approval of Newport Corporation. This document is provided for information only, and product specifications are subject to change without notice. Any change will be reflected in future publishings.

Table of Contents

1.0	Introduction	1
1.1	Purpose	1
1.2	Overview	1
<hr/>		
2.0	Command Interface.....	2
2.1	Constructor	2
2.2	Functions	2
2.2.1	General	2
	◆ CloseInstrument	2
	◆ GetDevices	2
	◆ OpenInstrument.....	2
	◆ WriteToInstrument	3
2.2.2	Commands.....	3
	◆ GP.....	3
	◆ ID_Get.....	3
	◆ ID_Set	4
	◆ IS_Get	4
	◆ IS_Set.....	4
	◆ IX_Get.....	5
	◆ IX_Set	5
	◆ IY_Get.....	5
	◆ IY_Set	6
	◆ LF_Get	6
	◆ LF_Set	6
	◆ OF_Get.....	7
	◆ OF_Set.....	7
	◆ PS_Get.....	8
	◆ PS_Set	8
	◆ PX_Get.....	8
	◆ PX_Set.....	9
	◆ PY_Get.....	9
	◆ PY_Set.....	9
	◆ PW_Get.....	10
	◆ PW_Set.....	10
	◆ RA	10
	◆ RC	11
	◆ RS.....	11
	◆ RS485.....	11

- ◆ SA_Get.....12
- ◆ SA_Set.....12
- ◆ TB_Get.....12
- ◆ TE.....13
- ◆ TS.....13
- ◆ VE.....13

3.0 Python Example..... 14

Service Form 17



Two-Axis Position & Power Sensing Device CONEX-PSD

1.0 Introduction

1.1 Purpose

The purpose of this document is to provide the method syntax of each command to communicate with the CONEX-PSD device.

1.2 Overview

The Command Interface is the wrapper class that maintains a list of CONEX-PSD instruments. It exposes methods to communicate with any CONEX-PSD device.

NOTE

Each function name is defined with the command code “AA”.

For each command function, refer to the CONEX-PSD programmer’s manual.

2.0 Command Interface

2.1 Constructor

ConexPSD()

The constructor is used to create an instance of the CONEX-PSD device.

2.2 Functions

2.2.1 General

◆ CloseInstrument

Syntax

int CloseInstrument()

return: 0 = successful or -1 = failure

Description

This function allows closing communication with the selected device. If the closing failed, the returned code is -1.

◆ GetDevices

Syntax

string[] GetDevices()

return: list of connected devices available to communicate

Description

This function returns the list of connected devices available to communicate.

◆ OpenInstrument

Syntax

int OpenInstrument(string strDeviceKey)

string strDeviceKey: device key

return: 0 = successful or -1 = failure

Description

This function allows opening communication with the selected device. If the opening failed, the returned code is -1.

◆ WriteToInstrument

Syntax

int WriteToInstrument(string command, ref string response, int stage)

command: Instrument command

response: Response of the command

stage: Instrument Stage

return:

Description

This Overridden function Queries or writes the command given by the user to the instrument.

2.2.2 Commands

◆ GP

Syntax

int GP(int controllerAddress, out double PositionX, out double PositionY, out double LaserPower, out string errstring)

controllerAddress: controllerAddress

PositionX: PositionX

PositionY: PositionY

LaserPower: LaserPower

errString: The failure reason

return: 0 in success and -1 on failure

Description

This function is used to process synchronous GP Get command which is used to Get X, Y positions and laser power level.

◆ ID_Get

Syntax

int ID_Get(int controllerAddress, out string SensorIdentifier, out string errstring)

controllerAddress: controllerAddress

SensorIdentifier: SensorIdentifier

errString: The failure reason

return: 0 in success and -1 on failure

Description

This function is used to process synchronous ID Get command which is used to get sensor identifier.

◆ **ID_Set**

Syntax

int ID_Set(int controllerAddress, string SensorIdentifier, out string errstring)

controllerAddress: controllerAddress

SensorIdentifier: SensorIdentifier

errString: The failure reason

return: 0 in success and -1 on failure

Description

This function is used to process synchronous ID Set command which is used to Set sensor identifier.

◆ **IS_Get**

Syntax

int IS_Get(int controllerAddress, out double Offset, out string errstring)

controllerAddress: controllerAddress

Offset: Offset

errString: The failure reason

return: 0 in success and -1 on failure

Description

This function is used to process synchronous IS Get command which is used to Get offset on ADC input SUM.

◆ **IS_Set**

Syntax

int IS_Set(int controllerAddress, double Offset, out string errstring)

controllerAddress: controllerAddress

Offset: Offset

errString: The failure reason

return: 0 in success and -1 on failure

Description

This function is used to process synchronous IS Set command which is used to Set offset on ADC input SUM.

◆ IX_Get**Syntax**

int IX_Get(int controllerAddress, out double OffsetADC1, out string errstring)

controllerAddress: controllerAddress

OffsetADC1: OffsetADC1

errString: The failure reason

return: 0 in success and -1 on failure

Description

This function is used to process synchronous IX Get command which is used to get offset on ADC input X.

◆ IX_Set**Syntax**

int IX_Set(int controllerAddress, double OffsetADC1, out string errstring)

controllerAddress: controllerAddress

OffsetADC1: OffsetADC1

errString: The failure reason

return: 0 in success and -1 on failure

Description

This function is used to process synchronous IX Set command which is used to set offset on ADC input X.

◆ IY_Get**Syntax**

int IY_Get(int controllerAddress, out double OffsetADC2, out string errstring)

controllerAddress: controllerAddress

OffsetADC2: OffsetADC2

errString: The failure reason

return: 0 in success and -1 on failure

Description

This function is used to process synchronous IY Get command which is used to get offset on ADC input Y.

◆ **IY_Set**

Syntax

int IY_Set(int controllerAddress, double OffsetADC2, out string errstring)

controllerAddress: controllerAddress

OffsetADC2: OffsetADC2

errString: The failure reason

return: 0 in success and -1 on failure

Description

This function is used to process synchronous IY Set command which is used to set offset on ADC input Y.

◆ **LF_Get**

Syntax

int LF_Get(int controllerAddress, out double Frequency, out string errstring)

controllerAddress: controllerAddress

Frequency: Frequency

errString: The failure reason

return: 0 in success and -1 on failure

Description

This function is used to process synchronous LF Get command which is used to get low pass filter frequency.

◆ **LF_Set**

Syntax

int LF_Set(int controllerAddress, double Frequency, out string errstring)

controllerAddress: controllerAddress

Frequency: Frequency

errString: The failure reason

return: 0 in success and -1 on failure

Description

This function is used to process synchronous LF Set command which is used to Set low pass filter frequency.

◆ OF_Get

NOTE

This command is not used with the CONEX-PSD9.

Syntax

int OF_Get(int controllerAddress, out double Offset1, out double Offset2, out double Offset3, out double Offset4, out string errstring)

controllerAddress: controllerAddress

Offset1: Offset #1

Offset2: Offset #2

Offset3: Offset #3

Offset4: Offset #4

errString: The failure reason

return: 0 in success and -1 on failure

Description

This function is used to process synchronous OF Get command which is used to get offsets. Refer to the CONEX-PSD Controller's manual to get the command description.

◆ OF_Set

NOTE

This command is not used with the CONEX-PSD9.

Syntax

int OF_Set(int controllerAddress, double Offset1, double Offset2, double Offset3, double Offset4, out string errstring)

controllerAddress: Controller's address

Offset1: Offset #1

Offset2: Offset #2

Offset3: Offset #3

Offset4: Offset #4

errString: The failure reason

return: 0 in success and -1 on failure

Description

This function is used to process synchronous OF Set command which is used to set offsets. Refer to the CONEX-PSD Controller's manual to get the command description.

◆ PS_Get**Syntax**

int PS_Get(int controllerAddress, out double Gain, out string errstring)

controllerAddress: controllerAddress

Gain: Gain

errString: The failure reason

return: 0 in success and -1 on failure

Description

This function is used to process synchronous PS Get command which is used to Get gain on ADC input SUM.

◆ PS_Set**Syntax**

int PS_Set(int controllerAddress, double Gain, out string errstring)

controllerAddress: controllerAddress

Gain: Gain

errString: The failure reason

return: 0 in success and -1 on failure

Description

This function is used to process synchronous PS Set command which is used to Set gain on ADC input SUM.

◆ PX_Get**Syntax**

int PX_Get(int controllerAddress, out double GainADC1, out string errstring)

controllerAddress: controllerAddress

GainADC1: GainADC1

errString: The failure reason

return: 0 in success and -1 on failure

Description

This function is used to process synchronous PX Get command which is used to Get gain on ADC input X.

◆ PX_Set**Syntax**

int PX_Set(int controllerAddress, double GainADC1, out string errstring)

controllerAddress: controllerAddress

GainADC1: GainADC1

errString: The failure reason

return: 0 in success and -1 on failure

Description

This function is used to process synchronous PX Set command which is used to Set gain on ADC input X.

◆ PY_Get**Syntax**

int PY_Get(int controllerAddress, out double GainADC2, out string errstring)

controllerAddress: controllerAddress

GainADC2: GainADC2

errString: The failure reason

return: 0 in success and -1 on failure

Description

This function is used to process synchronous PY Get command which is used to Get gain on ADC input Y.

◆ PY_Set**Syntax**

int PY_Set(int controllerAddress, double GainADC2, out string errstring)

controllerAddress: controllerAddress

GainADC2: GainADC2

errString: The failure reason

return: 0 in success and -1 on failure

Description

This function is used to process synchronous PY Set command which is used to Set gain on ADC input Y.

◆ PW_Get**Syntax**

int PW_Get(int controllerAddress, out int ConfigurationState, out string errstring)

controllerAddress: controllerAddress

ConfigurationState: ConfigurationState

errString: The failure reason

return: 0 in success and -1 on failure

Description

This function is used to process synchronous PW Get command which is used to Enter/Leave CONFIGURATION state.

◆ PW_Set**Syntax**

int PW_Set(int controllerAddress, int ConfigurationState, out string errstring)

controllerAddress: controllerAddress

ConfigurationState : ConfigurationState

errString: The failure reason

return: 0 in success and -1 on failure

Description

This function is used to process synchronous PW Set command which is used to Enter/Leave CONFIGURATION state.

◆ RA**Syntax**

int RA(int controllerAddress, out double RawAnalogInput1, out double RawAnalogInput2, out string errstring)

controllerAddress: controllerAddress

RawAnalogInput1: RawAnalogInput1

RawAnalogInput2: RawAnalogInput2

errString: The failure reason

return: 0 in success and -1 on failure

Description

This function is used to process synchronous RA Get command which is used to get raw analog input values.

◆ **RC**

Syntax

int RC(int controllerAddress, out double CorrectedAnalogInput1, out double CorrectedAnalogInput2, out string errstring)

controllerAddress: controllerAddress

CorrectedAnalogInput1: CorrectedAnalogInput1

CorrectedAnalogInput2: CorrectedAnalogInput2

errString: The failure reason

return: 0 in success and -1 on failure

Description

This function is used to process synchronous RC Get command which is used to Get corrected analog input values.

◆ **RS**

Syntax

int RS(int controllerAddress, out string errstring)

controllerAddress: controllerAddress

errString: The failure reason

return: 0 in success and -1 on failure

Description

This function is used to process synchronous RS Set command which is used to Reset controller.

◆ **RS485**

Syntax

int RS485(int controllerAddress, out string errstring)

controllerAddress: controllerAddress

errString: The failure reason

return: 0 in success and -1 on failure

Description

This function is used to process synchronous RS485 Set command which is used to Reset controller's address to 1.

◆ SA_Get**Syntax**

int SA_Get(int controllerAddress, out int Address, out string errstring)

controllerAddress: controllerAddress

Address: Address

errString: The failure reason

return: 0 in success and -1 on failure

Description

This function is used to process synchronous SA Get command which is used to get controller's RS-485 address.

◆ SA_Set**Syntax**

int SA_Set(int controllerAddress, int Address, out string errstring)

controllerAddress: controllerAddress

Address : Address

errString: The failure reason

return: 0 in success and -1 on failure

Description

This function is used to process synchronous SA Set command which is used to Set controller's RS-485 address.

◆ TB_Get**Syntax**

int TB_Get(int controllerAddress, string inErrorCode, string outErrorCode, out string errstring)

controllerAddress: controllerAddress

inErrorCode: Input error code (optional)

outErrorCode: Output error description

errString: The failure reason

return: 0 in success and -1 on failure

Description

This function is used to process synchronous TB Get command which is used to Get command error string.

◆ **TE****Syntax**

int TE(int controllerAddress, out string LastCommandError, out string errstring)

controllerAddress: controllerAddress

LastCommandError: LastCommandError

errString: The failure reason

return: 0 in success and -1 on failure

Description

This function is used to process synchronous TE Get command which is used to Get last command error.

◆ **TS****Syntax**

int TS(int controllerAddress, out string ErrorCode, out string StatusCode, out string errstring)

controllerAddress: controllerAddress

ErrorCode: ErrorCode

StatusCode: StatusCode

errString: The failure reason

return: 0 in success and -1 on failure

Description

This function is used to process synchronous TS Get command which is used to Get positioner error and controller state.

◆ **VE****Syntax**

int VE(int controllerAddress, out string ControllerVersion, out string errstring)

controllerAddress: controllerAddress

ControllerVersion: ControllerVersion

errString: The failure reason

return: 0 in success and -1 on failure

Description

This function is used to process synchronous VE Get command which is used to Get controller revision information.

3.0 Python Example

```

=====
#Initialization Start
#The script within Initialization Start and Initialization End is needed for properly
#initializing Command Interface for Conex-PSD instrument.
#The user should copy this code as is and specify correct paths here.
import sys

#Command Interface DLL can be found here.
print "Adding location of Newport.CONEXPSD.CommandInterface.dll to sys.path"
sys.path.append(r'C:\Program Files (x86)\Newport\MotionControl\CONEX-PSD\Bin')

# The CLR module provide functions for interacting with the underlying
# .NET runtime
import clr
# Add reference to assembly and import names from namespace
clr.AddReferenceToFile("Newport.CONEXPSD.CommandInterface.dll")
from CommandInterface import *

import System
=====
# Instrument Initialization
# The key should have double slashes since
# (one of them is escape character)
instrument="CONEX-PSD (A6T7NSPR)"
print 'Instrument Key=>', instrument

# create a device instance
PSD = ConexPSD()

componentID = PSD. OpenInstrument(instrument);
print 'componentID=>', componentID

# Get analog output #1 value
result, X, Y, LaserPower, errString = PSD.GP_Get(1)
if result == 0 :
    print 'X =>', X
    print 'Y =>', Y
    print 'Laser power =>', LaserPower
else:
    print 'Error=>',errString

```

```
# Get controller revision information
result, response, errString = PSD.VE(1)
if result == 0 :
    print 'controller revision=>', response
else:
    print 'Error=>',errString

# Get last command error
result, response, errString = PSD.TE(1)
if result == 0 :
    print 'Last command error =>', response
else:
    print 'Error=>',errString

# Unregister device
PSD. CloseInstrument();
```




Visit Newport Online at:
www.newport.com

North America & Asia

Newport Corporation
1791 Deere Ave.
Irvine, CA 92606, USA

Sales

Tel.: (800) 222-6440
e-mail: sales@newport.com

Technical Support

Tel.: (800) 222-6440
e-mail: tech@newport.com

Service, RMAs & Returns

Tel.: (800) 222-6440
e-mail: service@newport.com

Europe

MICRO-CONTROLE Spectra-Physics S.A.S
9, rue du Bois Sauvage
91055 Évry CEDEX
France

Sales

Tel.: +33 (0)1.60.91.68.68
e-mail: france@newport.com

Technical Support

e-mail: tech_europe@newport.com

Service & Returns

Tel.: +33 (0)2.38.40.51.55

