

843-R / 843-R-USB Field Upgrade Instructions


Reason for upgrade:

- 1.16
 - Added Japanese interface.
- 1.24
 - Added Russian and Chinese interfaces
 - Add support for 843-R-USB


Instructions

Field Upgrade of the 843-R / 843-R-USB firmware is a very simple process

1. After downloading the 843-R Field Upgrade Tool and copying it to your local drive, double-click on the file to run.
2. The 843-R Field Upgrade Tool expects .NET Framework 4 to be installed on the computer.

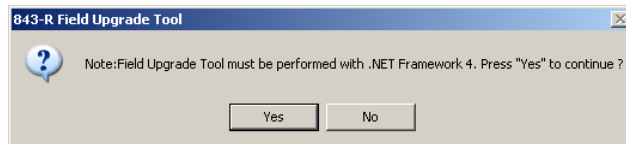

3. If not already installed, download .NET Framework 4 from Microsoft's web site at <http://www.microsoft.com/en-us/download/details.aspx?id=17851>
4. While the Field Upgrade Tool is self-extracting itself, the following will be displayed on the screen

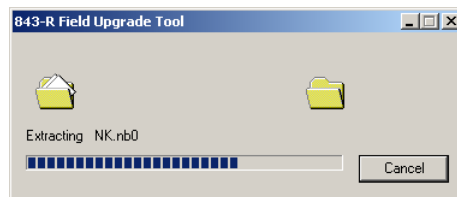

5. Follow the on-screen instructions to set up the 843-R / 843-R-USB for the firmware version upgrade.

