



Newport

Newport Power Meter

Firmware Upgrade Directions

Version 1.3

Revision Date: September 14, 2007

To Upgrade Firmware:

1. Extract all the files from PM1918-XXX.zip file. “XXX” denotes the version of firmware files present in the zip file. Password for the zip file is “newport”.
2. Put new firmware files (PM1918App.exe and XMLFile1.xml) in the root directory of a Win CE compatible USB flash drive (thumb drive).
3. Plug the USB flash drive into the USB port located on the front panel of a power meter, and turn it ON.
4. After initialization, the new firmware files will be detected and the power meter will ask if you want to download the files.
5. Press the “Yes” softkey to start the upgrade process.
6. The power meter will instruct you to restart once the upgrade is successful.
7. Restart the power meter by turning it OFF and back ON.
8. The power meter will restart running the new firmware.

Files in Firmware that can be updated:

PM1918App.exe

Main Firmware file

XMLFile1.xml

XML Screen Layout definitions