



# CONEX-CC

Single-Axis DC Motion  
with Controller/Driver



 **Newport®**

**LabVIEW Drivers**

**V2.0.x**

©2018 by Newport Corporation, Irvine, CA. All rights reserved.

Original instructions.

No part of this document may be reproduced or copied without the prior written approval of Newport Corporation. This document is provided for information only, and product specifications are subject to change without notice. Any change will be reflected in future publishings.

# Table of Contents

Preface.....	<a href="#">vi</a>
Confidentiality & Proprietary Rights .....	<a href="#">vi</a>
<b>1.0 CONEX-CC LabVIEW Drivers.....</b>	<b>1</b>
<b>2.0 VI's Libraries.....</b>	<b>4</b>
2.1 CONEX-ALL General Menu – Communication VI's .....	4
2.1.1   _Create Instrument Channel.vi .....	5
2.1.2   _Close Communications.vi.....	5
2.1.3   _Parse Instrument Channel.vi.....	6
2.1.4   Connect Newport Instrument.vi.....	6
2.1.5   FindInstrument.vi .....	7
2.1.6   NP_Logging.vi .....	7
2.2 CONEX-ALL Enter-Leave CONFIGURATION State v2.vi .....	8
2.3 CONEX-ALL Get Command Error String v2.vi .....	8
2.4 CONEX-ALL Get Controller Version v2.vi.....	8
2.5 CONEX-ALL Get Controller's address v2.vi .....	8
2.6 CONEX-ALL Get Identifier v2.vi.....	8
2.7 CONEX-ALL Get Last Command Error v2.vi.....	9
2.8 CONEX-ALL Get Positioner Error And Controller State v2.vi .....	9
2.9 CONEX-ALL Reset Controller v2.vi .....	9
2.10 CONEX-ALL Reset Controller's Address To 1 v2.vi .....	9
2.11 CONEX-ALL Set Controller's address v2.vi .....	9
2.12 CONEX-ALL Set Identifier v2.vi .....	9
2.13 Examples .....	10
2.14 CONEX Configuration.....	10
2.15 Motion Setup.....	10
2.16 Motion.....	11
2.17 Example .....	11
<b>3.0 VI'S Description .....</b>	<b>12</b>
3.1 CONEX-CC Absolute Move v2.vi .....	12
3.2 CONEX-CC Configure Simultaneous Started Move v2.vi.....	12
3.3 CONEX-CC Enter-Leave CONFIGURATION State v2.vi.....	13
3.4 CONEX-CC Enter-Leave DISABLE State v2.vi .....	14
3.5 CONEX-CC Execute Simultaneous Started Move v2.vi.....	14
3.6 CONEX-CC Get Acceleration v2.vi.....	14
3.7 CONEX-CC Get Backlash Compensation v2.vi.....	14

3.8	CONEX-CC Get Command Error String v2.vi .....	15
3.9	CONEX-CC Get Control Loop State v2.vi.....	15
3.10	CONEX-CC Get Controller Version v2.vi.....	15
3.11	CONEX-CC Get Controller's RS485 address v2.vi .....	16
3.12	CONEX-CC Get Current Configuration Parameters v2.vi.....	16
3.13	CONEX-CC Get Current Position v2.vi .....	16
3.14	CONEX-CC Get Derivative Gain v2.vi .....	16
3.15	CONEX-CC Get Driver Voltage v2.vi.....	17
3.16	CONEX-CC Get Encoder Increment Value v2.vi.....	17
3.17	CONEX-CC Get Following Error Limit v2.vi .....	17
3.18	CONEX-CC Get Friction Compensation v2.vi .....	18
3.19	CONEX-CC Get Home Search Time-out v2.vi .....	18
3.20	CONEX-CC Get Home Search Type v2.vi.....	18
3.21	CONEX-CC Get Home Search Velocity v2.vi .....	19
3.22	CONEX-CC Get Hysteresis Compensation v2.vi .....	19
3.23	CONEX-CC Get Integral Gain v2.vi .....	19
3.24	CONEX-CC Get Jerk Time v2.vi.....	20
3.25	CONEX-CC Get Last Command Error v2.vi.....	20
3.26	CONEX-CC Get Low Pass Filter for Kd v2.vi .....	20
3.27	CONEX-CC Get Motion Time for a Relative Move v2.vi.....	21
3.28	CONEX-CC Get Motor's Current Limits v2.vi .....	21
3.29	CONEX-CC Get Negative Software Limit v2.vi .....	22
3.30	CONEX-CC Get Positioner Error And Controller State v2.vi .....	22
3.31	CONEX-CC Get Positive Software Limit v2.vi.....	22
3.32	CONEX-CC Get Proportional Gain v2.vi .....	22
3.33	CONEX-CC Get SetPoint Position v2.vi .....	22
3.34	CONEX-CC Get Stage Identifier v2.vi .....	23
3.35	CONEX-CC Get Velocity Feed Forward v2.vi.....	23
3.36	CONEX-CC Get Velocity v2.vi.....	23
3.37	CONEX-CC Home search v2.vi .....	23
3.38	CONEX-CC Relative Move v2.vi.....	24
3.39	CONEX-CC Reset Controller v2.vi .....	24
3.40	CONEX-CC Reset Controller's Address To 1 v2.vi.....	24
3.41	CONEX-CC Set Acceleration v2.vi.....	24
3.42	CONEX-CC Set Backlash Compensation v2.vi.....	25
3.43	CONEX-CC Set Control Loop State v2.vi.....	25
3.44	CONEX-CC Set Controller's RS-485 address v2.vi.....	25
3.45	CONEX-CC Set Derivative Gain v2.vi.....	26
3.46	CONEX-CC Set Driver Voltage v2.vi .....	26
3.47	CONEX-CC Set Encoder Increment Value v2.vi .....	26
3.48	CONEX-CC Set Following Error Limit v2.vi .....	26
3.49	CONEX-CC Set Friction Compensation v2.vi .....	27
3.50	CONEX-CC Set Home Search Time-out v2.vi .....	27
3.51	CONEX-CC Set Home Search Type v2.vi.....	27

---

3.52	CONEX-CC Set Home Search Velocity v2.vi .....	27
3.53	CONEX-CC Set Hysteresis Compensation v2.vi.....	28
3.54	CONEX-CC Set Integral Gain v2.vi .....	28
3.55	CONEX-CC Set Jerk Time v2.vi .....	28
3.56	CONEX-CC Set Low Pass Filter for Kd v2.vi .....	28
3.57	CONEX-CC Set Motor's Current Limits v2.vi.....	29
3.58	CONEX-CC Set Negative Software Limit v2.vi.....	29
3.59	CONEX-CC Set Positive Software Limit v2.vi .....	29
3.60	CONEX-CC Set Proportional Gain v2.vi.....	30
3.61	CONEX-CC Set Stage Identifier v2.vi.....	30
3.62	CONEX-CC Set Velocity Feed Forward v2.vi .....	30
3.63	CONEX-CC Set Velocity v2.vi.....	30
3.64	CONEX-CC Stop Motion v2.vi .....	31
	<b>Service Form .....</b>	<b>33</b>

---

# Preface

## Confidentiality & Proprietary Rights

### **Reservation of Title**

The Newport Programs and all materials furnished or produced in connection with them ("Related Materials") contain trade secrets of Newport and are for use only in the manner expressly permitted. Newport claims and reserves all rights and benefits afforded under law in the Programs provided by Newport Corporation.

Newport shall retain full ownership of Intellectual Property Rights in and to all development, process, align or assembly technologies developed and other derivative work that may be developed by Newport. Customer shall not challenge, or cause any third party to challenge, the rights of Newport.

### **Preservation of Secrecy and Confidentiality and Restrictions to Access**

Customer shall protect the Newport Programs and Related Materials as trade secrets of Newport, and shall devote its best efforts to ensure that all its personnel protect the Newport Programs as trade secrets of Newport Corporation. Customer shall not at any time disclose Newport's trade secrets to any other person, firm, organization, or employee that does not need (consistent with Customer's right of use hereunder) to obtain access to the Newport Programs and Related Materials. These restrictions shall not apply to information (1) generally known to the public or obtainable from public sources; (2) readily apparent from the keyboard operations, visual display, or output reports of the Programs; (3) previously in the possession of Customer or subsequently developed or acquired without reliance on the Newport Programs; or (4) approved by Newport for release without restriction.

## Sales, Tech Support & Service

### **North America & Asia**

Newport Corporation  
1791 Deere Ave.  
Irvine, CA 92606, USA

#### **Sales**

Tel.: (800) 222-6440  
e-mail: sales@newport.com

#### **Technical Support**

Tel.: (800) 222-6440  
e-mail: tech@newport.com

#### **Service, RMAs & Returns**

Tel.: (800) 222-6440  
e-mail: service@newport.com

### **Europe**

MICRO-CONTROLE Spectra-Physics S.A.S  
9, rue du Bois Sauvage  
91055 Evry Cedex  
France

#### **Sales France**

Tel.: +33 (0)1.60.91.68.68  
e-mail: france@newport.com

#### **Technical Support**

e-mail: tech\_europe@newport.com

#### **Service & Returns**

Tel.: +33 (0)2.38.40.51.55

First printing 2010.

**Copyright 2010-2013** by Newport Corporation, Irvine, CA. All rights reserved. No part of this manual may be reproduced or copied without the prior written approval of Newport Corporation. This manual is provided for information only, and product specifications are subject to change without notice. Any change will be reflected in future printings.



# Single-Axis DC Motor Controller/Driver CONEX-CC


## 1.0 CONEX-CC LabVIEW Drivers

The CONEX-CC LabVIEW drivers have been developed under LABVIEW 2010

### NOTE

You must use at least the [2010](#) of LabVIEW.

Copy the directory CONEX-CC Controller\_Drivers under the directory **user.lib** (\CONEX\CONEX-CC) of LabVIEW 20xx.

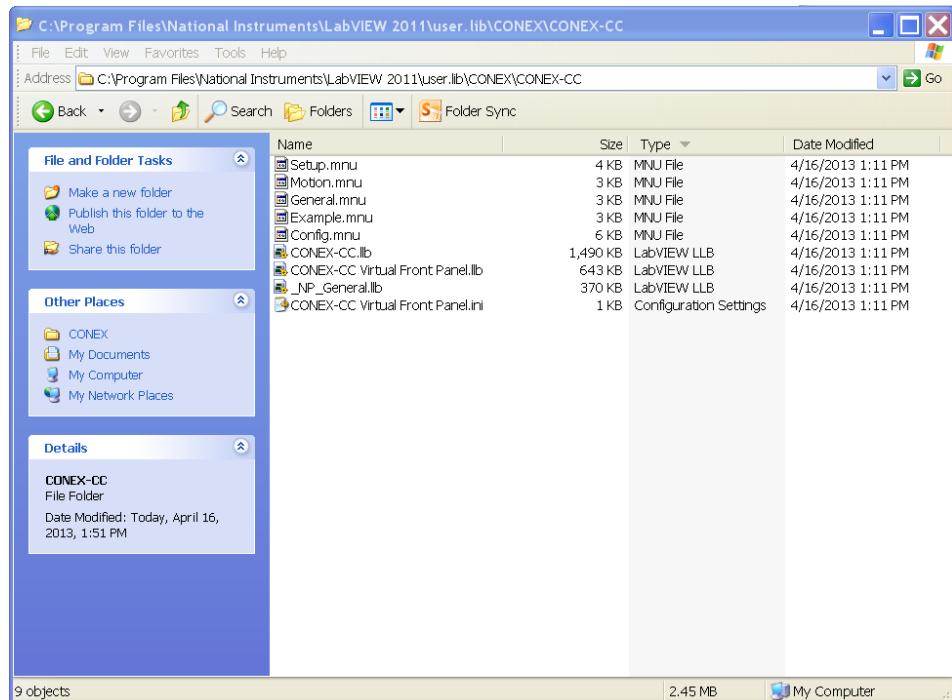

This directory contains documented VIs, menu to access the different VIs and controls defined to use the CONEX-CC, and the different menus where the VIs will be in LabVIEW:

**Config:** VIs to change configuration parameters (those used for setup after reset or end of configuration mode).

**Conex-ALL:** VIs to change configuration that all CONEX devices have and general communication Vis.

**Mot. Setup:** VIs to change working parameters (those lost when switching off the controller).

- Motion:** VIs to move or stop the positioner.  
**Example:** CONEX-CC Virtual Front panel VIs.

You select both **CONEX** menu and CONEX-CC Controller-Drivers sub-menu from User Libraries:



When you activate the Help window, you will see the description of each of the VIs.  
Click on a menu then select a VI. Place it and connect it.

You must use the connection VIs to setup connection. The Connect CONEX-CC device.vi will find the CONEX on the USB ports and setup connection for you, so that you can just connect it to the first subvi. These VIs are in the CONEX-ALL (General) menu.

Communication settings:

COM port – (Use the samples to find it or look under Device Manager)  
Baud Rate – 921600 (USB serial speed)

In each VI, there is a Communication Cluster that contains the following elements:

**VISA resource name in:** VISA resource name is passed to low level VI's

**Device name in:** Readable description of device

**Controller address in:** Channel number

#### Note on Controller Address:

(Important for other devices with multiple RS485 connections, used to match command syntax of these similar RS485 instruments, however, for USB connection only one channel is addressed per USB cable, so all addresses can be set to 1, regardless of number of connected devices.)

**error in:** The **error in** cluster can accept error information wired from VIs previously called. Use this information to decide if any functionality should be bypassed in the event of errors from other VIs.

The pop-up option **Explain Error** (or Explain Warning) gives more information about the error displayed.

**status:** The **code** input identifies the error or warning.

The pop-up option **Explain Error** (or Explain Warning) gives more information about the error displayed.

**code:** The code input identifies the error or warning.

The pop-up option **Explain Error** (or Explain Warning) gives more information about the error displayed.



**source:** The **source** string describes the origin of the error or warning.

The pop-up option **Explain Error** (or Explain Warning) gives more information about the error displayed.


## 2.0 VI's Libraries

### 2.1

#### CONEX-ALL General Menu – Communication VI's

The Communication VIs at the top are low-level sub-VIs that talk to the device for you. The lower VIs with white background are configuration Vis common to all the CONEX family of devices.

The “Connect Newport Instrument.vi” will setup a connection and build a Connection Cluster that is all you need to pass to the other CONEX-CC specific function sub-VIs.

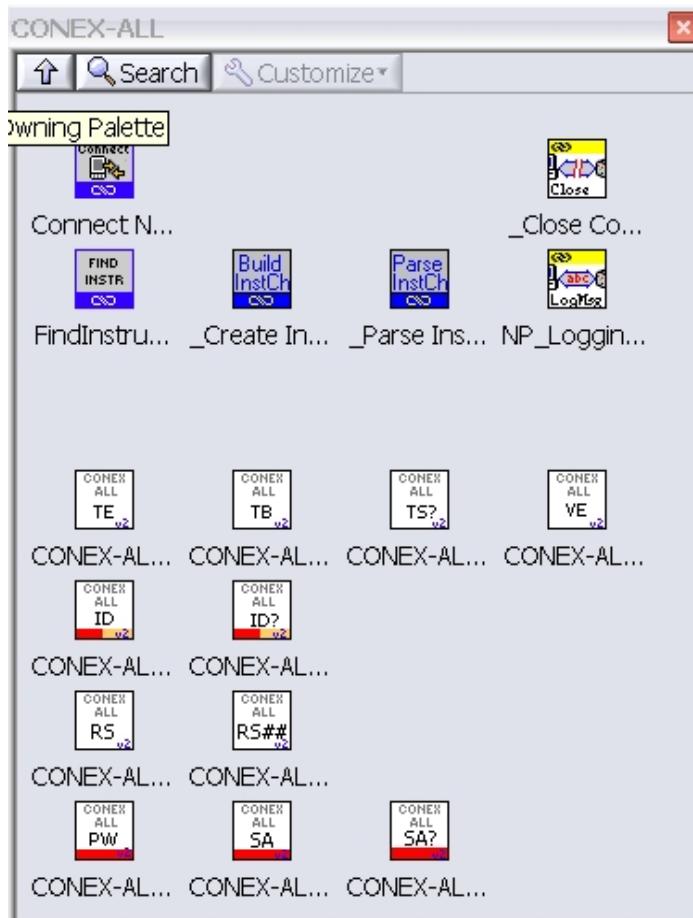


### 2.1.1 \_Create Instrument Channel.vi

Builds the instrument channel handle (string) for an instrument connected over GPIB or RS-232. The default output String for a USB is RS-232 at a Baud Rate of 921600.

This string should be passed in and out of the library VIs to control the specified instrument. Different handles should be used for controlling multiple instruments.

This will also initialize the RS-232 port to the speed specified; which must be done manually if not using this VI.

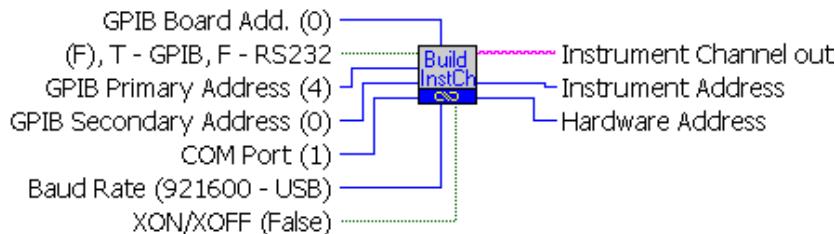

**[TFI]** (F), T - GPIB, F - RS232

**[U8]** GPIB Primary Address (4)

**[U16]** COM Port (1)

**[U8]** GPIB Secondary Address (0)

**[U8]** GPIB Board Add. (0)

**[U32]** Baud Rate (921600 - USB)


**[TFI]** XON/XOFF (False)

**[abc]** Instrument Channel out

**[I16]** Instrument Address

**[I32]** Hardware Address

### 2.1.2 \_Close Communications.vi

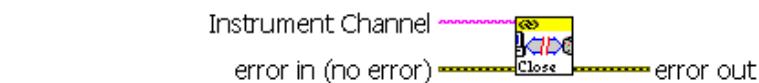


**[abc]** error in (no error)

**[abc]** Instrument Channel

**[abc]** Instrument Channel

### 2.1.3 Parse Instrument Channel.vi

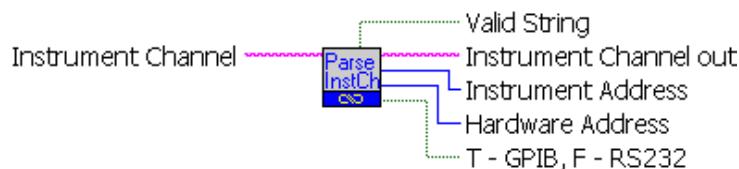

**Instrument Channel in**

**Instrument Channel out**

**Instrument Address**

**Hardware Address**

**T - GPIB, F - RS232**


**Valid String**

### 2.1.4 Connect Newport Instrument.vi

Connect Newport Device

Get list of instrument, filter (if set) and allow for selection of device to talk to.

This is generic selection of the instrument to connected to, so look in system settings or on device to verify it is the correct port.

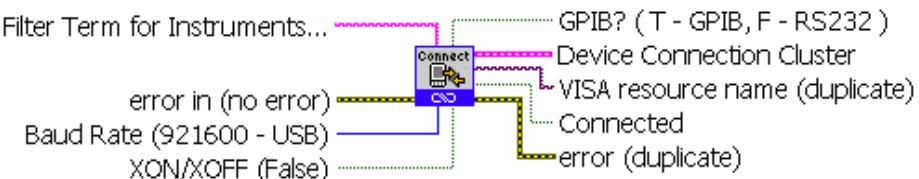

**Instrument Filter Term for Instruments (" - none)**  
String used to verify expected device is found.

**error in (no error)**

**Baud Rate (921600 - USB)**

**XON/XOFF (False)**


**Connected**  
Connected = true when connection is successful

**VISA resource name (duplicate)**

**error (duplicate)**

**GPIB? ( T - GPIB, F - RS232 )**

### 2.1.5 FindInstrument.vi

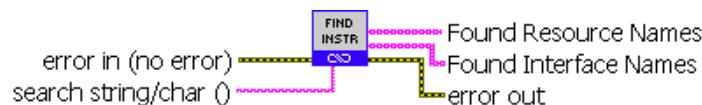

**search string/char ()**


**error in (no error)**

**error out**

**Found Resource Names**

**Found Interface Names**

### 2.1.6 NP\_Logging.vi

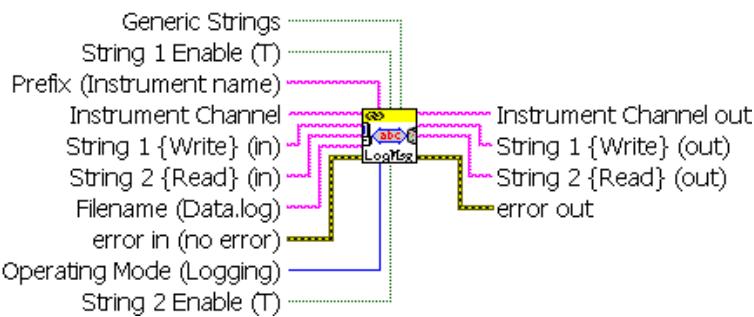

**error in (no error)**

**Operating Mode (Logging)**

**String 1 {Write} (in)**

**String 2 {Read} (in)**

**Filename (Data.log)**

**Generic Strings**

**String 1 Enable (T)**

**String 2 Enable (T)**

**Prefix (Instrument name)**

**Instrument Channel**

**error out**

**String 2 {Read} (out)**

**String 1 {Write} (out)**

**Instrument Channel out**

## 2.2 CONEX-ALL Enter-Leave CONFIGURATION State v2.vi




**[TF]** **Go to CONFIGURATION State**  
**Configuration State?**  
**T - In Configuration State**

## 2.3 CONEX-ALL Get Command Error String v2.vi



**[abc]** **Error code**  
**Error Code**  
**[abc]** **Error description**  
**Description of input error code.**

## 2.4 CONEX-ALL Get Controller Version v2.vi

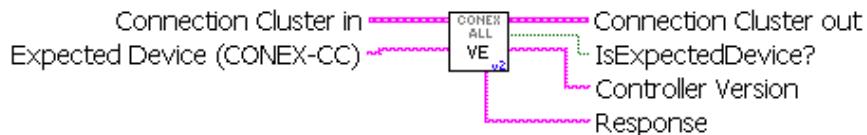

**[abc]** **Expected Device (CONEX-CC)**  
**[abc]** **Response**  
**Full Response**  
**[abc]** **Controller Version**  
**Controller Version information**  
**[TF]** **IsExpectedDevice?**  
**Is CONEX - CC?**

## 2.5 CONEX-ALL Get Controller's address v2.vi



**[abc]** **Controller's address**  
**Controllers address**

## 2.6 CONEX-ALL Get Identifier v2.vi




**[abc]** **Identifier**  
**Identification of attached hardware.**

## 2.7 CONEX-ALL Get Last Command Error v2.vi



**abc** **Error code**  
Last Command Error

## 2.8 CONEX-ALL Get Positioner Error And Controller State v2.vi

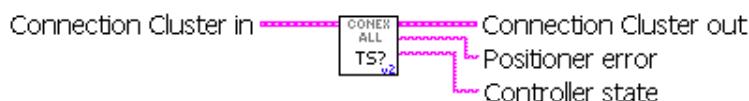

**abc** **Positioner error**  
Positioner Error

**abc** **Controller state**  
Controller State

## 2.9 CONEX-ALL Reset Controller v2.vi



Resets CONEX-CC

## 2.10 CONEX-ALL Reset Controller's Address To 1 v2.vi



## 2.11 CONEX-ALL Set Controller's address v2.vi



**132** **Controller's address**  
Controller's RS-485 address


## 2.12 CONEX-ALL Set Identifier v2.vi




**abc** **Stage Identifier**  
Stage Identifier

## 2.13 Examples


CONEX-Sample.vi shows how easy it is to find, connect and get version:

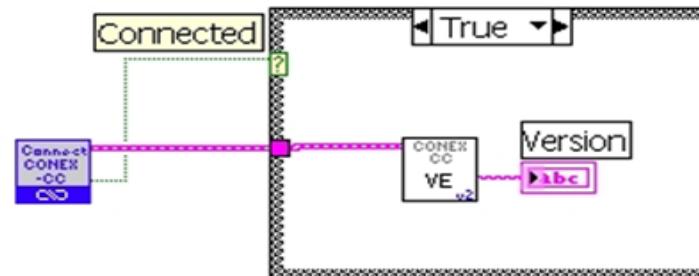


## 2.14 CONEX Configuration

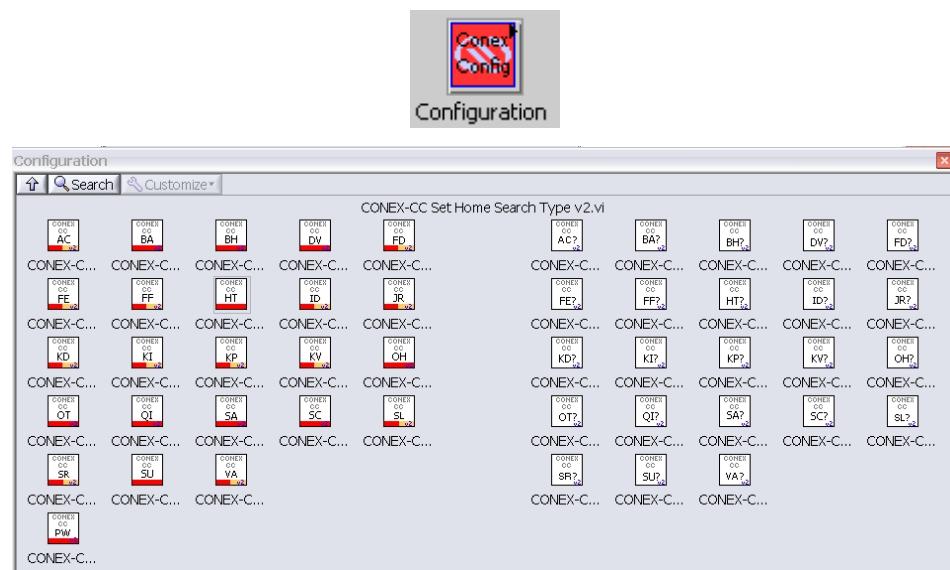


## 2.15 Motion Setup

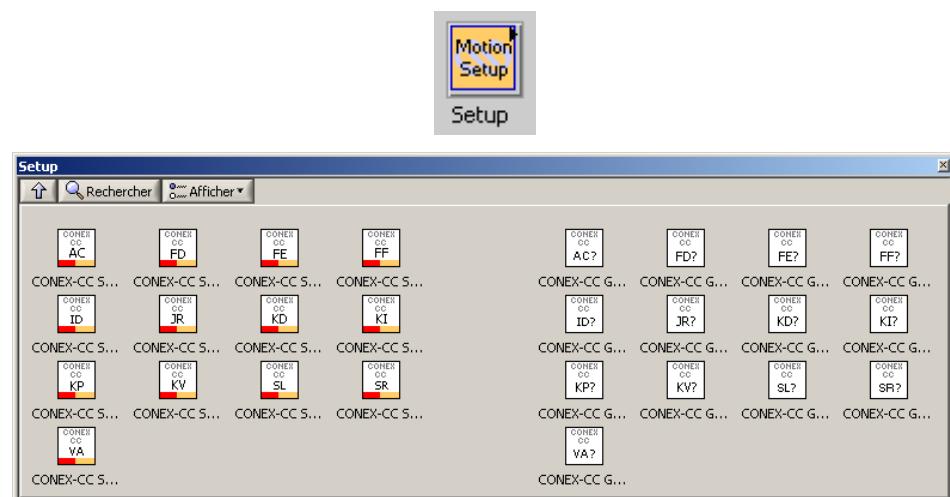


## 2.16 Motion

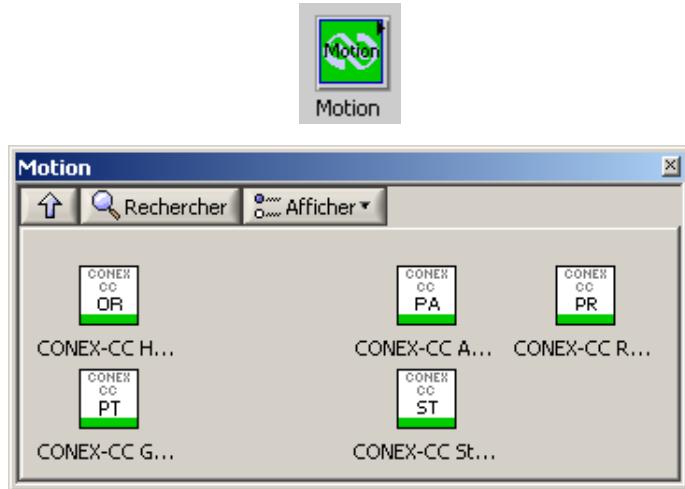


## 2.17 Example

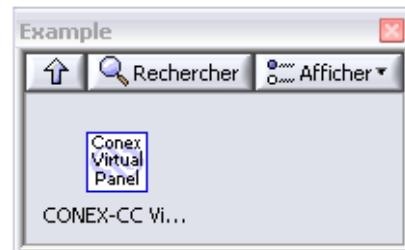

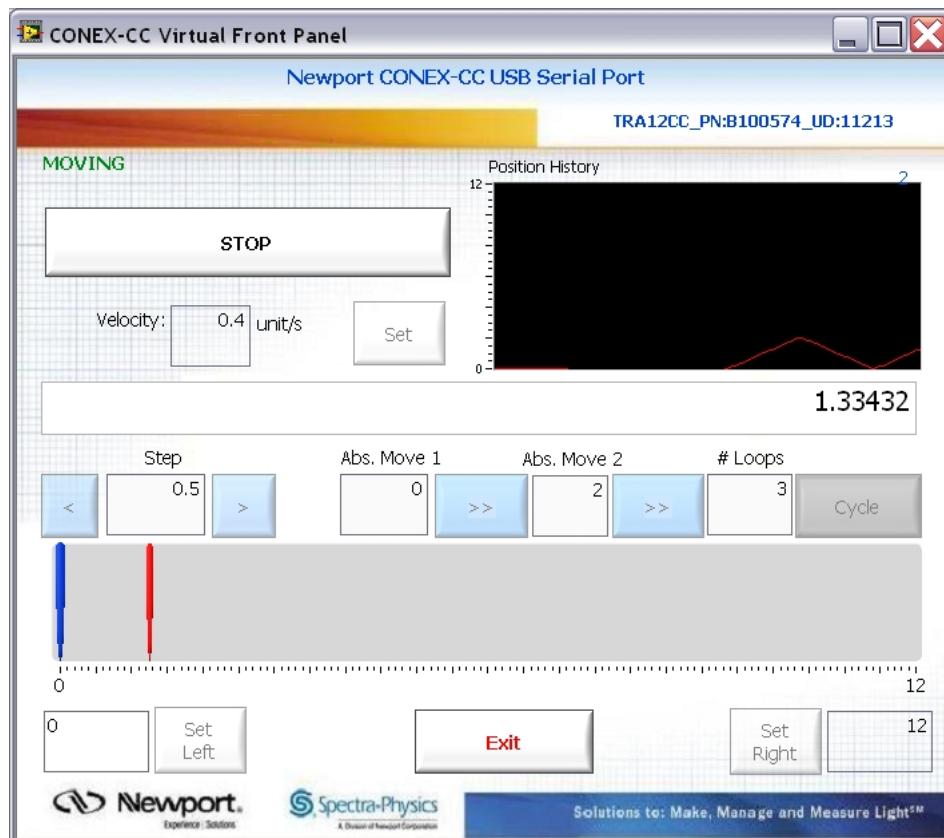

## 3.0 VI'S Description

### 3.1 CONEX-CC Absolute Move v2.vi

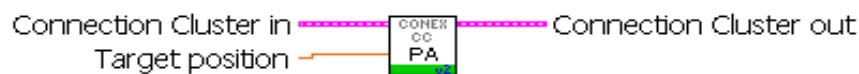

Set Absolute Position Move to start.

**Target position**  
Absolute Position target

### 3.2 CONEX-CC Configure Simultaneous Started Move v2.vi




Set Simultaneous Started Move.

**Target position**

Simultaneous Move Position target

### 3.3 CONEX-CC Enter-Leave CONFIGURATION State v2.vi

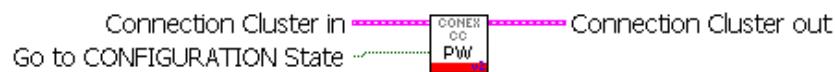

Set Configuration State to Enabled or Disabled.

T - Enabled - Enter Configuration state.


F - Leave Configuration state.

**Go to CONFIGURATION State**

Configuration State?

T - In Configuration State

### 3.4 CONEX-CC Enter-Leave DISABLE State v2.vi

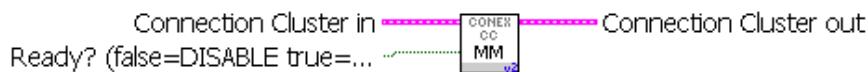

Enable Ready State or Disabled State

Generally used after connecting to allow for moves to start, as device starts in Disabled state.



**Ready? (false=DISABLE true=READY)**

Ready? flag  
true - Ready to move  
false – Disabled

### 3.5 CONEX-CC Execute Simultaneous Started Move v2.vi



**SE** - Execute simultaneous started move

### 3.6 CONEX-CC Get Acceleration v2.vi




**AC?** - Get acceleration



**Acceleration**

Acceleration value

### 3.7 CONEX-CC Get Backlash Compensation v2.vi

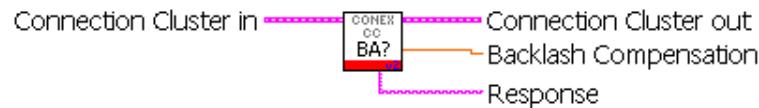

**BA?** - Get backlash compensation



**Response**


Full Backlash Compensation Response string



**Backlash Compensation**

Backlash Compensation

### 3.8 CONEX-CC Get Command Error String v2.vi

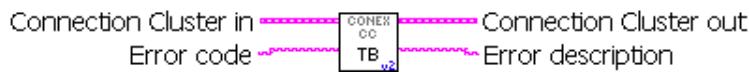


Get description of error from error code.

**abc** **Error code**  
Error Code

**abc** **Error description**  
Description of input error code.

### 3.9 CONEX-CC Get Control Loop State v2.vi

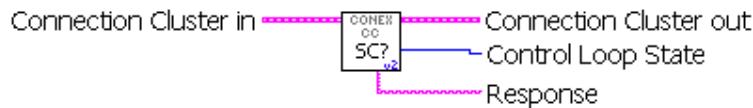


**SC?** - Get control loop state

**abc** **Response**  
Contol Loop State full Response

**I32** **Control Loop State**  
Contol Loop State

### 3.10 CONEX-CC Get Controller Version v2.vi

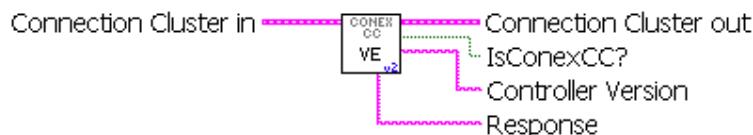

**VE** - Get controller revision information

**abc** **Controller Version**  
Controller Version information

**TF** **IsConexCC?**  
Is CONEX - CC?

### 3.11 CONEX-CC Get Controller's RS485 address v2.vi



**SA?** - Set controller's RS-485 address


### 3.12 CONEX-CC Get Current Configuration Parameters v2.vi



**ZT** - Get current configuration parameters

**abc** **Configuration parameters**  
Configuration parameters for device

### 3.13 CONEX-CC Get Current Position v2.vi

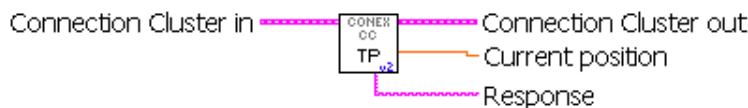


**TP?** - Get Current Position

**abc** **Response**  
Current Position Response

**DBL** **Current position**  
Current Position


### 3.14 CONEX-CC Get Derivative Gain v2.vi

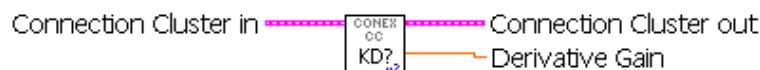

**KD?** - Get derivative gain

**DBL** **Derivative Gain**  
Derivative Gain

### 3.15 CONEX-CC Get Driver Voltage v2.vi

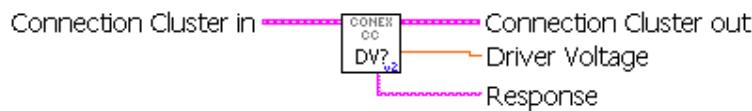


DV? - Get driver voltage

**Response**  
Driver Voltage

**Driver Voltage**  
Driver Voltage

### 3.16 CONEX-CC Get Encoder Increment Value v2.vi

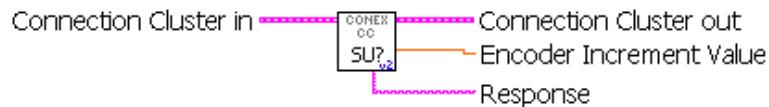


SU? - Get encoder increment value

**Response**  
Encoder Increment Value Response

**Encoder Increment Value**  
Encoder Increment Value

### 3.17 CONEX-CC Get Following Error Limit v2.vi

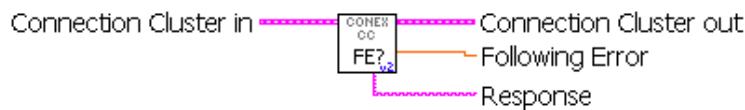


FE? - Get following error limit

**Response**  
Following Error Response

**Following Error**  
Following Error

### 3.18 CONEX-CC Get Friction Compensation v2.vi

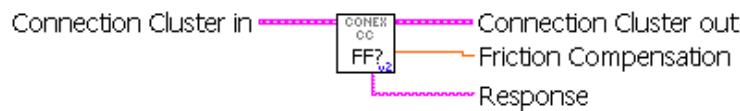


**FF?** - Get friction compensation

**abc** **Response**  
Friction Compensation Response

**DBL** **Friction Compensation**  
Friction Compensation

### 3.19 CONEX-CC Get Home Search Time-out v2.vi

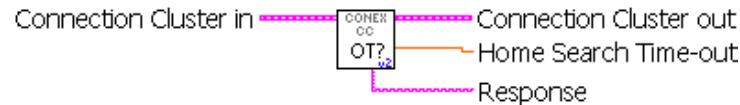


**OT?** - Get HOME search time-out

**abc** **Response**  
Home Search Time-out Response

**DBL** **Home Search Time-out**  
Home Search Time-out

### 3.20 CONEX-CC Get Home Search Type v2.vi

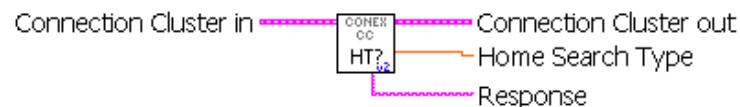


**HT?** - Get HOME search type

**abc** **Response**  
HOME search type Response

**DBL** **Home Search Type**  
HOME search type

### 3.21 CONEX-CC Get Home Search Velocity v2.vi

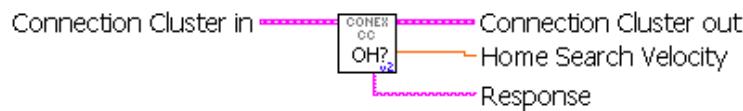


**OH?** - Get HOME search velocity

**abc** **Response**  
HOME search velocity Response

**DBL** **Home Search Velocity**  
HOME search velocity

### 3.22 CONEX-CC Get Hysteresis Compensation v2.vi

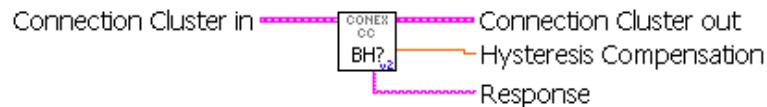

**BH?** - Get hysteresis compensation

**abc** **Response**  
Hysteresis compensation Response

**DBL** **Hysteresis Compensation**  
Hysteresis compensation


### 3.23 CONEX-CC Get Integral Gain v2.vi



**KI?** - Get integral gain

**DBL** **Integral Gain**  
Integral gain

### 3.24 CONEX-CC Get Jerk Time v2.vi

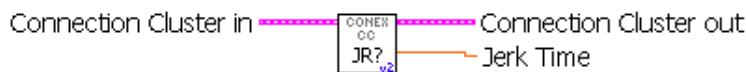

**JR?** - Get jerk time




### 3.25 CONEX-CC Get Last Command Error v2.vi





**TE?** - Get Error




### 3.26 CONEX-CC Get Low Pass Filter for Kd v2.vi

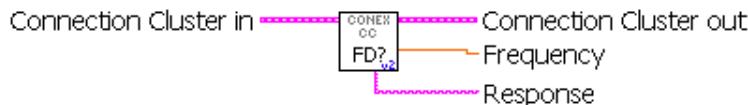

**FD?** - Get low pass filter cut off frequency for Kd

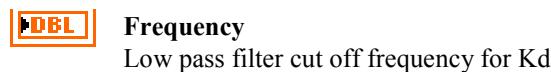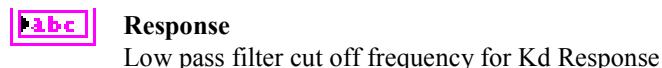

### 3.27 CONEX-CC Get Motion Time for a Relative Move v2.vi

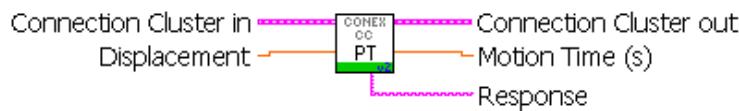

**PT** - Get motion time for a relative move

Input size of move to make, and it will output time for move.



#### Displacement

Displacement to find time to achieve



#### Response


Motion time for a relative move Response



#### Motion Time (s)

Motion time for a relative move

### 3.28 CONEX-CC Get Motor's Current Limits v2.vi

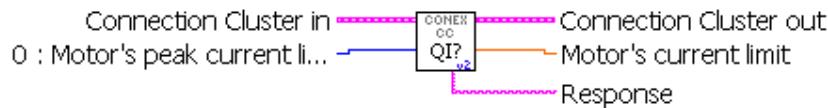

**QIx?** - Get motor's current limits

Motor peak current selection of limit to query:

0: Motor's peak current limit (default)

1: Motor's rms current limit

2: Motor's rms current averaging time



#### Motor Query Type (0 - Peak Current)



#### Response

Motor's current limits Response



#### Motor's current limit

Motor's current limits


### 3.29 CONEX-CC Get Negative Software Limit v2.vi



Get Negative Software Limit

**DBL** Motor Query Type (0 - Peak Current) Left limit  
Negative Software Limit

### 3.30 CONEX-CC Get Positioner Error And Controller State v2.vi

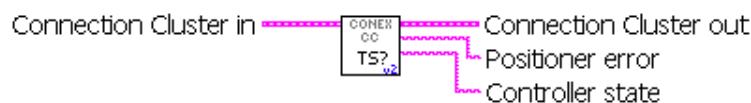


TS? - Get Positioner Error and Controller State

### 3.31 CONEX-CC Get Positive Software Limit v2.vi




SR? - Get Positive Software Limit

### 3.32 CONEX-CC Get Proportional Gain v2.vi

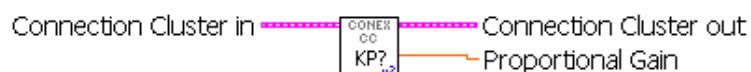

KP? - Get proportional gain


### 3.33 CONEX-CC Get SetPoint Position v2.vi

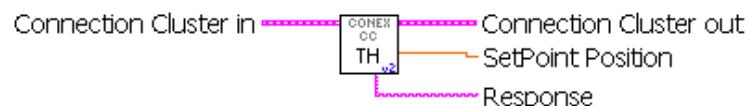


TH? - Get set-point position

**3.34 CONEX-CC Get Stage Identifier v2.vi**

**ID?** - Get Stage Identifier

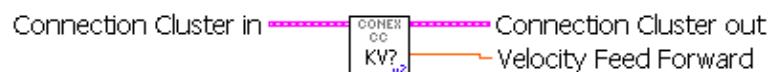**3.35 CONEX-CC Get Velocity Feed Forward v2.vi**


**KV?** - Get velocity feed forward

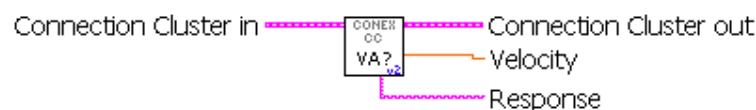**3.36 CONEX-CC Get Velocity v2.vi**


**VA?** - Get Velocity

**Response**  
Velocity Response


**Velocity**  
Velocity

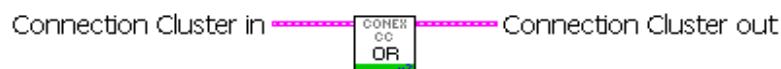**3.37 CONEX-CC Home search v2.vi**

Home stage

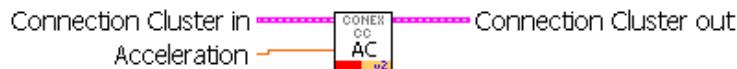**3.38 CONEX-CC Relative Move v2.vi****PR - Relative Move** **Relative displacement**  
Relative Move Displacement**3.39 CONEX-CC Reset Controller v2.vi****RS - Reset controller****3.40 CONEX-CC Reset Controller's Address To 1 v2.vi****RS## - Reset controller's address to 1****3.41 CONEX-CC Set Acceleration v2.vi****AC - Set acceleration** **Acceleration**  
Acceleration

### 3.42 CONEX-CC Set Backlash Compensation v2.vi

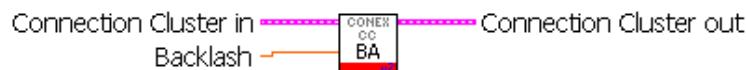

**BA** - Set backlash compensation

**[DBL]** **Backlash**  
Backlash compensation

### 3.43 CONEX-CC Set Control Loop State v2.vi



**SC** - Set control loop state

**[132]** **Control Loop State**  
Control loop state

### 3.44 CONEX-CC Set Controller's RS-485 address v2.vi



**SA** - Set controller's RS-485 address


---

#### NOTE

For CONEX over USB the RS-485 Address will always be 1, this command will not change the address. Each USB bus (cable) only talks to the attached device, not to any other device. Therefore, having the option of changing the address is not needed.


---

**[132]** **Controller's address**  
Controller's RS-485 address

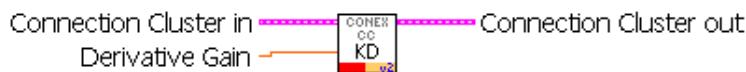**3.45 CONEX-CC Set Derivative Gain v2.vi**

**KD** - Set derivative gain

**DBL** **Derivative Gain**  
Derivative gain

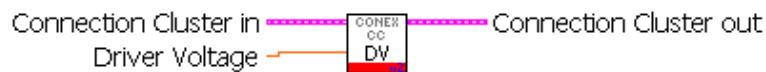**3.46 CONEX-CC Set Driver Voltage v2.vi**


**DV** - Set driver voltage

**DBL** **Driver Voltage**  
Driver voltage

**3.47 CONEX-CC Set Encoder Increment Value v2.vi**

**SU** - Set encoder increment value

**DBL** **Encoder increment value**  
Encoder increment value

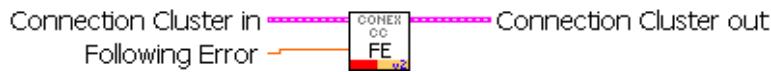**3.48 CONEX-CC Set Following Error Limit v2.vi**


**FE** - Set following error limit

**DBL** **Following Error**  
Following error limit


### 3.49 CONEX-CC Set Friction Compensation v2.vi




FF - Set friction compensation

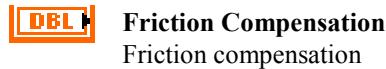


### 3.50 CONEX-CC Set Home Search Time-out v2.vi

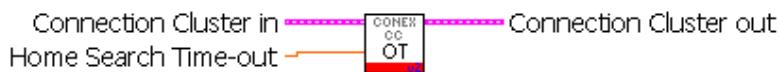

OT - Set Home Search Time-out

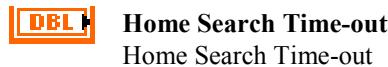


### 3.51 CONEX-CC Set Home Search Type v2.vi

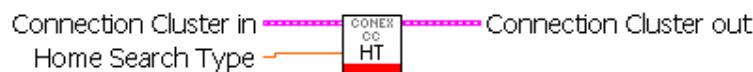


HT - Set HOME search type



### 3.52 CONEX-CC Set Home Search Velocity v2.vi

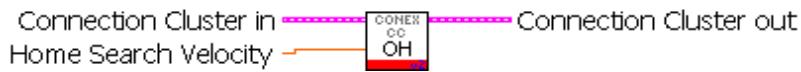

OH - Set HOME search velocity

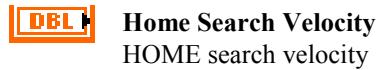


### 3.53 CONEX-CC Set Hysteresis Compensation v2.vi



**BH** - Set hysteresis compensation

**DBL** **Hysteresis**  
Hysteresis

### 3.54 CONEX-CC Set Integral Gain v2.vi

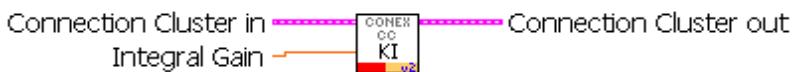

**KI** - Set integral gain

**DBL** **Integral Gain**  
Integral Gain


### 3.55 CONEX-CC Set Jerk Time v2.vi



**JR** - Set jerk time

**DBL** **Jerk Time**  
Jerk Time


### 3.56 CONEX-CC Set Low Pass Filter for Kd v2.vi

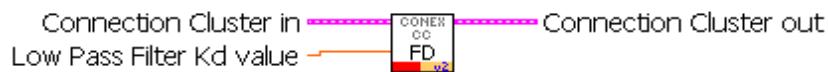

**FD** - Set Low Pass Filter for Kd

**DBL** **Low Pass Filter Kd value**  
Low Pass Filter Kd value

### 3.57 CONEX-CC Set Motor's Current Limits v2.vi

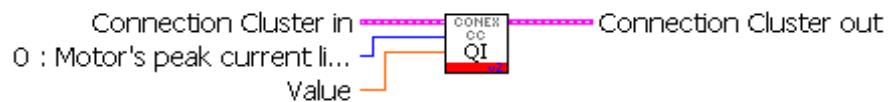

**QIx** - Set motor's current limits

Motor Limit Types:

0: Motor's peak current limit (default)

1: Motor's rms current limit

2: Motor's rms current averaging time

The Type of limit must be selected and the value to set.

#### NOTE

Right click on the Type and selecting {Create Constant} will create an easy selection for the type.


**[DBL]** **Value**

New Motor's Limit for Type selected

**[I32]** **Motor's Limit Type Select (0 - Peak current)**

Motor Limit Type select

### 3.58 CONEX-CC Set Negative Software Limit v2.vi

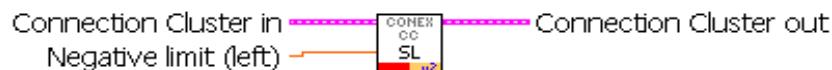

Set Negative Limit of travel

**[DBL]** **Negative limit (left)**

### 3.59 CONEX-CC Set Positive Software Limit v2.vi




**SR** - Set Positive limit (right)

**[DBL]** **Positive limit (right)**

Positive limit (right)


### 3.60 CONEX-CC Set Proportional Gain v2.vi

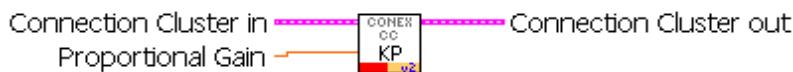

**KP** - Set proportional gain




### 3.61 CONEX-CC Set Stage Identifier v2.vi

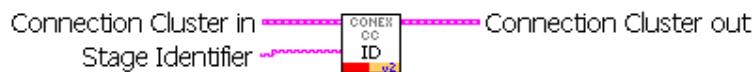


**ID** - Set Stage Identifier




### 3.62 CONEX-CC Set Velocity Feed Forward v2.vi

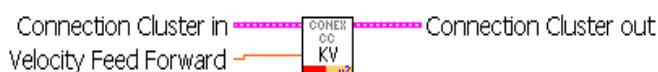

**KV** - Set velocity feed forward

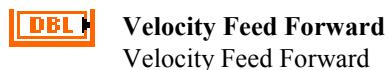

### 3.63 CONEX-CC Set Velocity v2.vi

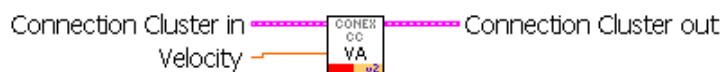

**VA** - Set Velocity



### 3.64 CONEX-CC Stop Motion v2.vi



**ST** - Stop Motion



## Service Form

Your Local Representative

Tel.: \_\_\_\_\_

Fax: \_\_\_\_\_

Name: \_\_\_\_\_

Return authorization #: \_\_\_\_\_

*(Please obtain prior to return of item)*

Company: \_\_\_\_\_

Date: \_\_\_\_\_

Address: \_\_\_\_\_

Phone Number: \_\_\_\_\_

Country: \_\_\_\_\_

Fax Number: \_\_\_\_\_

P.O. Number: \_\_\_\_\_

Serial #: \_\_\_\_\_

Item(s) Being Returned: \_\_\_\_\_

Model#: \_\_\_\_\_

Description: \_\_\_\_\_

Reasons of return of goods (please list any specific problems): \_\_\_\_\_

---

---

---

---

---

---

---

---

---

---

---

---

---

---

---

---

---

---

---

---

---



Visit Newport Online at:

[www.newport.com](http://www.newport.com)

**North America & Asia**

Newport Corporation  
1791 Deere Ave.  
Irvine, CA 92606, USA

**Sales**

Tel.: (800) 222-6440  
e-mail: [sales@newport.com](mailto:sales@newport.com)

**Technical Support**

Tel.: (800) 222-6440  
e-mail: [tech@newport.com](mailto:tech@newport.com)

**Service, RMAs & Returns**

Tel.: (800) 222-6440  
e-mail: [service@newport.com](mailto:service@newport.com)

**Europe**

MICRO-CONTROLE Spectra-Physics S.A.S  
9, rue du Bois Sauvage  
91055 Évry CEDEX  
France

**Sales**

Tel.: +33 (0)1.60.91.68.68  
e-mail: [france@newport.com](mailto:france@newport.com)

**Technical Support**

e-mail: [tech\\_europe@newport.com](mailto:tech_europe@newport.com)

**Service & Returns**

Tel.: +33 (0)2.38.40.51.55



**Newport®**

**Ophir®**

**Spectra-Physics®**