



## FTP (File Transfer Protocol) Connection to the XPS with Internet Explorer (IE) Version 7 and Later


If using the ftp://XPS IP Address results in an error from IE indicating that the page could not be displayed:

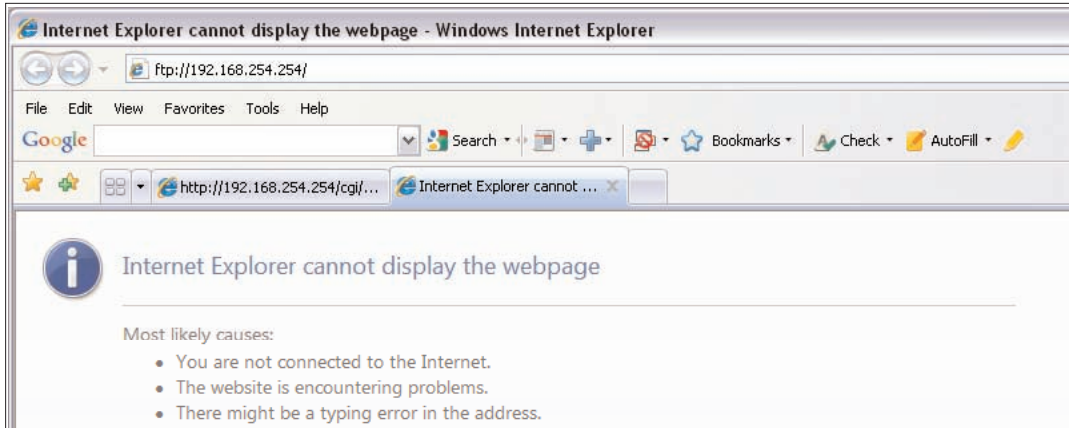

1. In the menu bar click on “Page” and select “Open FTP site in Windows Explorer”.

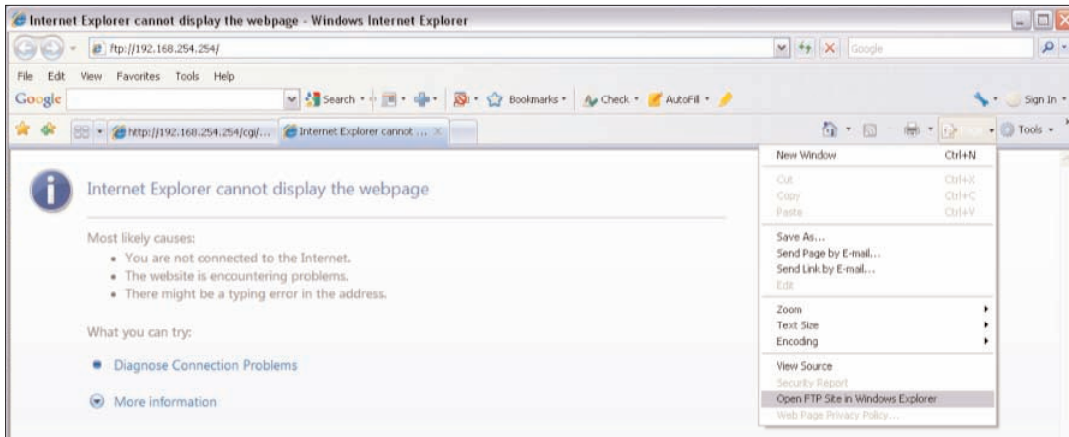

2. A new IE window will open and the following “FTP Folder Error” will display.


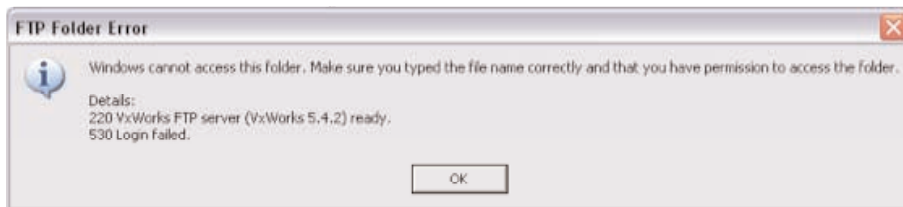


## FTP (File Transfer Protocol) Connection to the XPS with Internet Explorer (IE) Version 7 and Later

3. Click on OK and select "File" and then "Login As..." from the menu bar.



4. Login using your username and password.

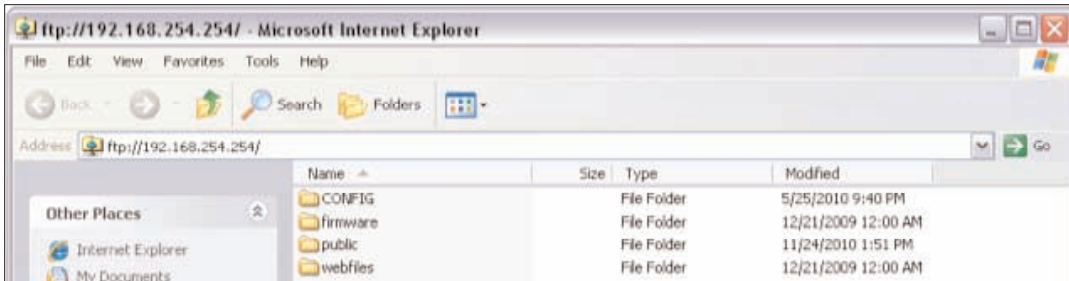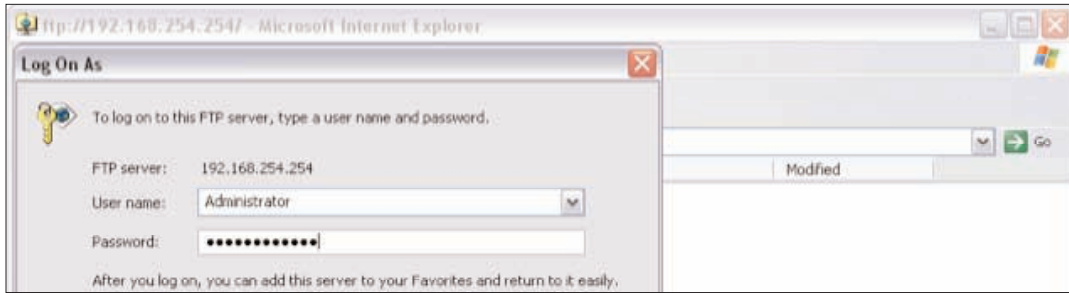

For more information please contact Newport Corporation Application Engineers at 800.222.6440.



**Newport Corporation, Global Headquarters**  
1791 Deere Avenue, Irvine, CA 92606, USA

[www.newport.com](http://www.newport.com)

**PHONE:** 1-800-222-6440 1-949-863-3144 **FAX:** 1-949-253-1680 **EMAIL:** [sales@newport.com](mailto:sales@newport.com)

Complete listings for all global office locations are available online at [www.newport.com/contact](http://www.newport.com/contact)

	PHONE
Belgium	+32-(0)0800-11 257
China	+86-10-6267-0065
France	+33-(0)1-60-91-68-68
Japan	+81-3-3794-5511
Taiwan	+886 -(0)2-2508-4977

EMAIL
<a href="mailto:belgium@newport.com">belgium@newport.com</a>
<a href="mailto:china@newport.com">china@newport.com</a>
<a href="mailto:france@newport.com">france@newport.com</a>
<a href="mailto:spectra-physics@splasers.co.jp">spectra-physics@splasers.co.jp</a>
<a href="mailto:sales@newport.com.tw">sales@newport.com.tw</a>

	PHONE
Irvine, CA, USA	+1-800-222-6440
Netherlands	+31-(0)30 6592111
United Kingdom	+44-1235-432-710
Germany / Austria / Switzerland	+49-(0)6151-708-0

EMAIL
<a href="mailto:sales@newport.com">sales@newport.com</a>
<a href="mailto:netherlands@newport.com">netherlands@newport.com</a>
<a href="mailto:uk@newport.com">uk@newport.com</a>
<a href="mailto:germany@newport.com">germany@newport.com</a>