

APEX2

Oriel[®] Illumination Systems



APEX2-XE

APEX2-QTH

User's Manual

Oriel[®]
INSTRUMENTS
A Newport Corporation Brand

TABLE OF CONTENTS

1	GENERAL INFORMATION	5
1.1	SYMBOLS AND DEFINITIONS	5
1.2	GENERAL WARNINGS	6
1.3	ELECTRICAL HAZARDS	6
1.4	FIRE HAZARDS	7
1.5	LAMP HANDLING	7
2	INTRODUCTION	8
3	SYSTEM SETUP	13
3.1	ITEMS INCLUDED WITH SYSTEM	13
3.2	UNPACKING	13
3.3	CHOOSING A LOCATION	13
3.4	ASSEMBLY	14
3.5	LAMP INSTALLATION	17
3.6	ELECTRICAL CONNECTION	24
3.7	LAMP INTENSITY ADJUSTMENT (APEX2-QTH ONLY)	26
4	CONTROLLING THE APEX2 FILTER WHEEL	27
4.1	MANUAL CONTROL	28
4.2	DIRECT APEX2 COMPUTER CONTROL	28
4.3	MONOCHROMATOR UTILITY SOFTWARE CONTROL	29
4.4	MONOCHROMATOR COMMAND SET	31
4.5	MONOCHROMATOR HAND CONTROLLER	31
4.6	TRACQ BASIC CONTROL	32
5	APEX2 FEATURES	33
5.1	FILTER WHEEL OVERVIEW	33
5.2	SHUTTER	33
5.3	ADJUSTABLE COLLIMATING LENS	34
5.4	MATING FLANGE OPTIONS	35
6	FILTER INSTALLATION	36
7	UNIVERSAL FILTER WHEEL UTILITY SOFTWARE	38
7.1	SYSTEM REQUIREMENTS	38
7.2	SOFTWARE INSTALLATION PROCEDURE	39
7.3	USING THE FILTER WHEEL UTILITY SOFTWARE	43
7.4	COMMAND SET	45
8	ACCESSORIES	46
8.1	MONOCHROMATOR MOUNTING KITS	46
8.2	FIBER FOCUSING ASSEMBLIES	46
8.3	FOCUSING LENS ASSEMBLIES	47
8.4	BEAM TURNERS	47
8.5	UV SAFETY ACCESSORIES	48
9	MAINTENANCE	49
9.1	LAMP REPLACEMENT	49
9.2	CLEANING	49
9.3	LAMP CARE AND HANDLING	49
9.4	FUSE REPLACEMENT	49
10	SPECIFICATIONS	50
10.1	APEX2 ILLUMINATOR	50
10.2	UNIVERSAL FILTER WHEEL UTILITY SOFTWARE	51
11	EU DECLARATION OF CONFORMITY	52

12	WARRANTY AND SERVICE	53
12.1	CONTACTING ORIEL INSTRUMENTS	53
12.2	REQUEST FOR ASSISTANCE/ SERVICE.....	54
12.3	REPAIR SERVICE	54
12.4	NON-WARRANTY REPAIR	54
12.5	WARRANTY REPAIR	55
12.6	LOANER/ DEMO MATERIAL.....	56

LIST OF FIGURES

Figure 1:	Basic System Block Diagram.....	9
Figure 2:	Dimensional Diagram	10
Figure 3:	APEX2-XE Lamp Spectra	11
Figure 4:	APEX2-QTH Lamp Spectra	12
Figure 5:	How the APEX2 unit arrives in its packaging.....	14
Figure 6:	Front lens assembly and integrated light source.....	15
Figure 7:	Female flange of the integrated light source unit (left) and installed front lens assembly (right).....	16
Figure 8:	Installed front lens assembly with Collimating Lens Adjustment Ring set screw oriented vertically for easy accessibility	16
Figure 9:	Remove Top Cover Screws	17
Figure 10:	Remove Housing Cover.....	18
Figure 11:	Remove Lamp Cable Holder	18
Figure 12:	Move Lamp Holder Wires	19
Figure 13:	Unfasten Power Supply Cable	19
Figure 14:	Insert Lamp	20
Figure 15:	Lamp Ellipse Positioning Groove	20
Figure 16:	Secure Lamp.....	21
Figure 17:	Connect Lamp Wiring	21
Figure 18:	Re-attach Lamp Cable Holder.....	22
Figure 19:	Completed Lamp Installation	22
Figure 20:	Safety Interlock	23
Figure 21:	Cover Re-installed.....	23
Figure 22:	APEX2 Power Connection	25
Figure 23:	Lamp Intensity Adjustment Settings.....	26
Figure 24:	Filter Wheel Connectors	27
Figure 25:	Manual Control Configuration	28
Figure 26:	Direct Computer Control Configuration.....	28
Figure 27:	Monochromator Control Configuration	29
Figure 28:	MS257 Monochromator Connections	30
Figure 29:	Cornerstone USB Monochromator Connections.....	30
Figure 30:	APEX2 Configured with Hand Controller	31
Figure 31:	Filter Selection Table in TracQ Basic	32
Figure 32:	Shutter Knob	33
Figure 33:	Collimating Lens Adjustment	34
Figure 34:	Do Not Adjust Relay Lens.....	34
Figure 35:	Oriel 1.5-Inch Series Flange	35
Figure 36:	Removing the Front Panel of the Filter Wheel.....	36
Figure 37:	Filter Wrench Removing Retaining Ring	37
Figure 38:	Installed Filters	37
Figure 39:	Setup Application	39
Figure 40:	Destination Directory Selection.....	39
Figure 41:	Accept License Agreements	40

Figure 42: Start Installation	40
Figure 43: Finish Installation	41
Figure 44: Restart Computer.....	41
Figure 45: Start Menu and Desktop Changes.....	42
Figure 46: Safely removing the USB stick	42
Figure 47: Connect to Filter Wheel	43
Figure 48: Device ID Query.....	43
Figure 49: Select Filter	44
Figure 50: Query Filter Position	44
Figure 51: Fiber Focusing Assembly	46
Figure 52: Model 6195 Optic Holder	47
Figure 53: Beam Turning Reflector Holder	47
Figure 54: Protective Gloves	48
Figure 55: 49125 and 49126 UV Safety Glasses	48
Figure 56: Fuse Compartment	49

1 GENERAL INFORMATION

Thank you for your purchase of this APEX2 illumination system from Oriel Instruments.

Please carefully read the following important safety precautions prior to unpacking and operating this equipment. In addition, please read the complete User's Manual for additional important notes and cautionary statements regarding the use and operation of the system.

1.1 SYMBOLS AND DEFINITIONS

	WARNING Situation has the potential to cause bodily harm or death.
	CAUTION Situation has the potential to cause damage to property or equipment.
	ELECTRICAL SHOCK HAZARD Hazard arising from dangerous voltage. Any mishandling could result in irreparable damage to the equipment, and personal injury or death.
	EUROPEAN UNION CE MARK The presence of the CE Mark on Newport Corporation equipment means that it has been designed, tested and certified as complying with all applicable European Union (CE) regulations and recommendations.
NOTE:	Additional important information the user or operator should consider.

Please read all instructions that were provided prior to operation of the system.

If there are any questions, please contact Oriel Instruments or the representative through whom the system was purchased.

1.2 GENERAL WARNINGS

- Read all warnings and operating instructions for this system prior to setup and use.
- Do not use this equipment in or near water.
- To prevent damage to the equipment, read the instructions in the equipment manual for proper input voltage.
- This equipment is grounded through the grounding conductor of the power cord.
- Route the power cord and other cables so they are not likely to be damaged.
- Disconnect power before cleaning the equipment.
- Do not use liquid or aerosol cleaners; use only a dry lint-free cloth.
- Lock out all electrical power sources before servicing the equipment.
- To avoid explosion, do not operate this equipment in an explosive atmosphere.
- Qualified service personnel should perform safety checks after any service.
- If this equipment is used in a manner not specified in this manual, the protection provided by this equipment may be impaired.
- To prevent damage to equipment when replacing fuses, locate and correct the problem that caused the fuse to blow before re-applying power.
- Do not block ventilation openings.
- Do not position this product in such a manner that would make it difficult to disconnect the power cord.
- Use only the specified replacement lamp.
- This product should only be powered as described in the manual.
- Do not remove the cover for normal usage.

1.3 ELECTRICAL HAZARDS

- Make all connections to or from the power supply with the power off.
- Do not use the power supply without its cover in place. Lethal voltages are present inside.

1.4 FIRE HAZARDS

Lamps are extremely hot during operation, and for several minutes after being shut off. Keep flammable objects away from the lamp and lamp housing.

Newport Research (fan-cooled) Housings are equipped with a condenser lens. The re-focused output of this lens can cause ignition of flammable targets including but not limited to walls, certain chemicals.

To avoid fire hazard, use only the specified fuses with the correct type number, voltage and current ratings as referenced in the appropriate locations in the service instructions or on the equipment. Only qualified service personnel should replace fuses.

1.5 LAMP HANDLING

Read all information and warnings provided with lamp.

The Xenon arc lamp used in the APEX2-XE is filled with rare gas at high pressure, presenting the danger of lamp explosion due to mechanical failure. This is particularly true when the lamp is operating, as the internal pressure can reach tens of atmospheres. Thermal strains can also cause the lamp to explode under certain conditions.

Never touch a lamp's glass envelope or the reflector's inner surface with bare fingers or other contaminants. Skin oil or other substances can burn into the lamp envelope during operation and negatively affect the lamp's performance and lifetime.

Always wear appropriate powder free gloves and impact-resistant goggles when handling any lamp. Avoid any mechanical strain during handling. Do not operate the lamp without all housing panels in place.

Lamps become very hot after only a few minutes of operation (up to 150°C) and remain quite hot for at least 10 to 15 minutes after being turned off.

2 INTRODUCTION

The APEX2 illumination systems are designed to be compact, economical, fully integrated solutions suitable for a wide variety of applications. The APEX2 is a plug and play light source producing a collimated 1.3 inch (33 mm) diameter output beam. Simply install the lamp, connect the power cord and switch on the power – then the system is ready for use.

The systems are available with either a Xenon or Quartz Tungsten Halogen (QTH) lamp. Lamp replacement is easy – no alignment is needed. The lamp operates with a built-in power supply, so no current or power presets or adjustments are required. The APEX2 is designed with an ellipsoidal reflector, capturing additional light to increase optical throughput.

The APEX2 is designed to be quick and easy to set up. Mounting feet help with handling and leveling the illuminator. The mounting feet are removable, so the illuminator may be permanently mounted to an optical table if desired. The APEX2 comes preassembled on a plate with mounting holes for securing the illuminator to an inch or metric optical table / breadboard.

The light source is coupled to a filter wheel, where up to six neutral density or ordering sorting filters may be installed. A large, easy to read digital display indicates which filter is in the light path. Filters are sold separately, as their specifications are application-dependent.

The filters may be selected using a variety of methods, including manual control and software. For those who wish to create their own programs, low level commands may be sent to the system via USB connection from a personal computer (PC).

The APEX2 includes a manually operated shutter. It is important that the lamp is warmed up and stabilized when using the system. When the lamp is not needed, it can remain on (and therefore warmed up) by closing off the light path.

NOTE: Always allow the lamp to warm up for at least 30 minutes prior to beginning work.

The output flange of the APEX2 is compatible with the 1.5-inch Oriel flanging system. The collimated output may be coupled to a wide variety of items such as a beam turner, focusing optics, fiber optic cable adapter or detector.

Compatible instruments:

- Cornerstone 130B Monochromator and its predecessor Cornerstone 130 Monochromator
- Cornerstone 260 Monochromator
- MS260i Spectrograph
- MS257 Monochromator
- MS257 Spectrograph

NOTE:

In this user's manual, the term "monochromator" is used throughout the text. The APEX2 may be operated with both monochromators and spectrographs.

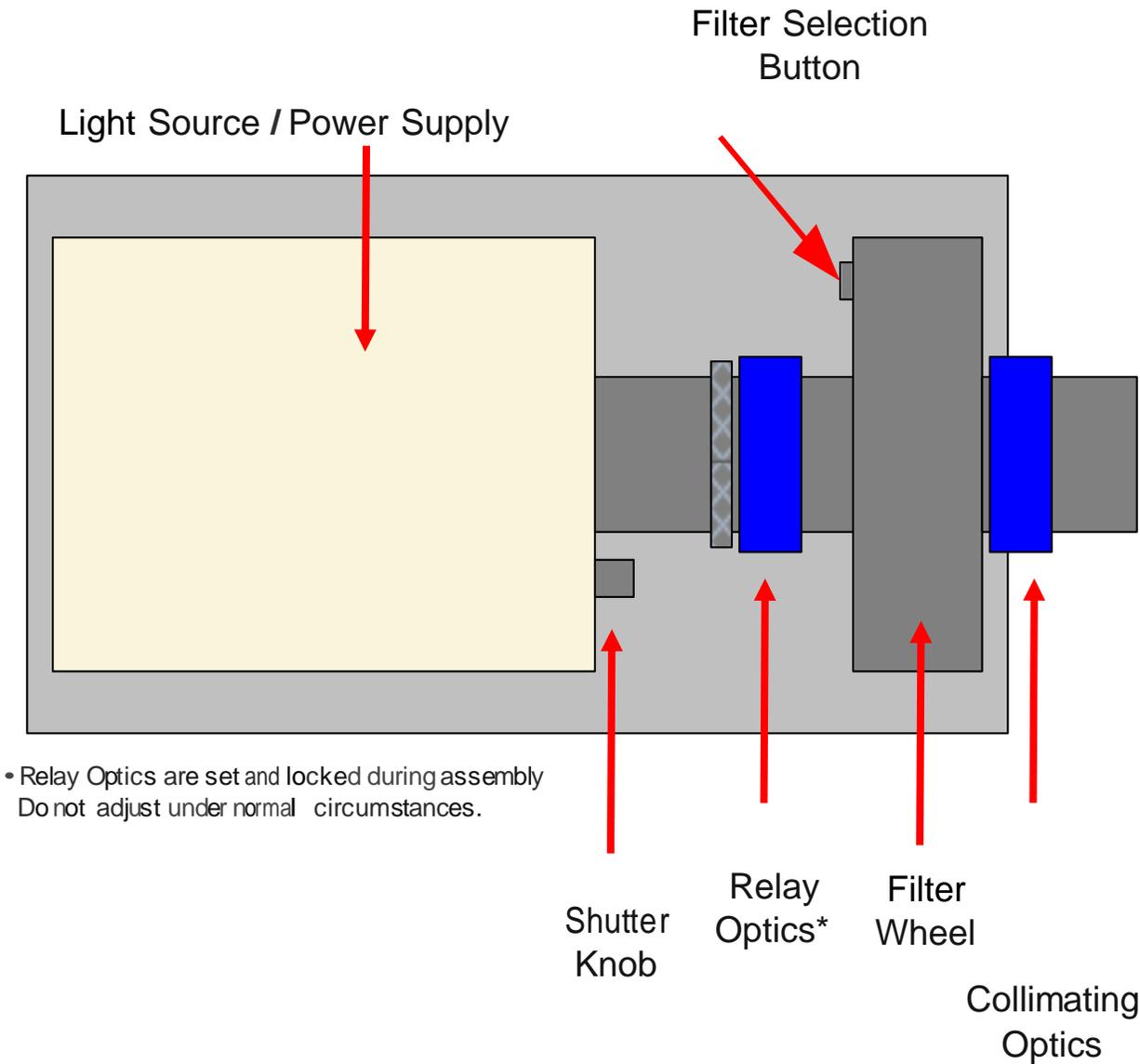


Figure 1: Basic System Block Diagram

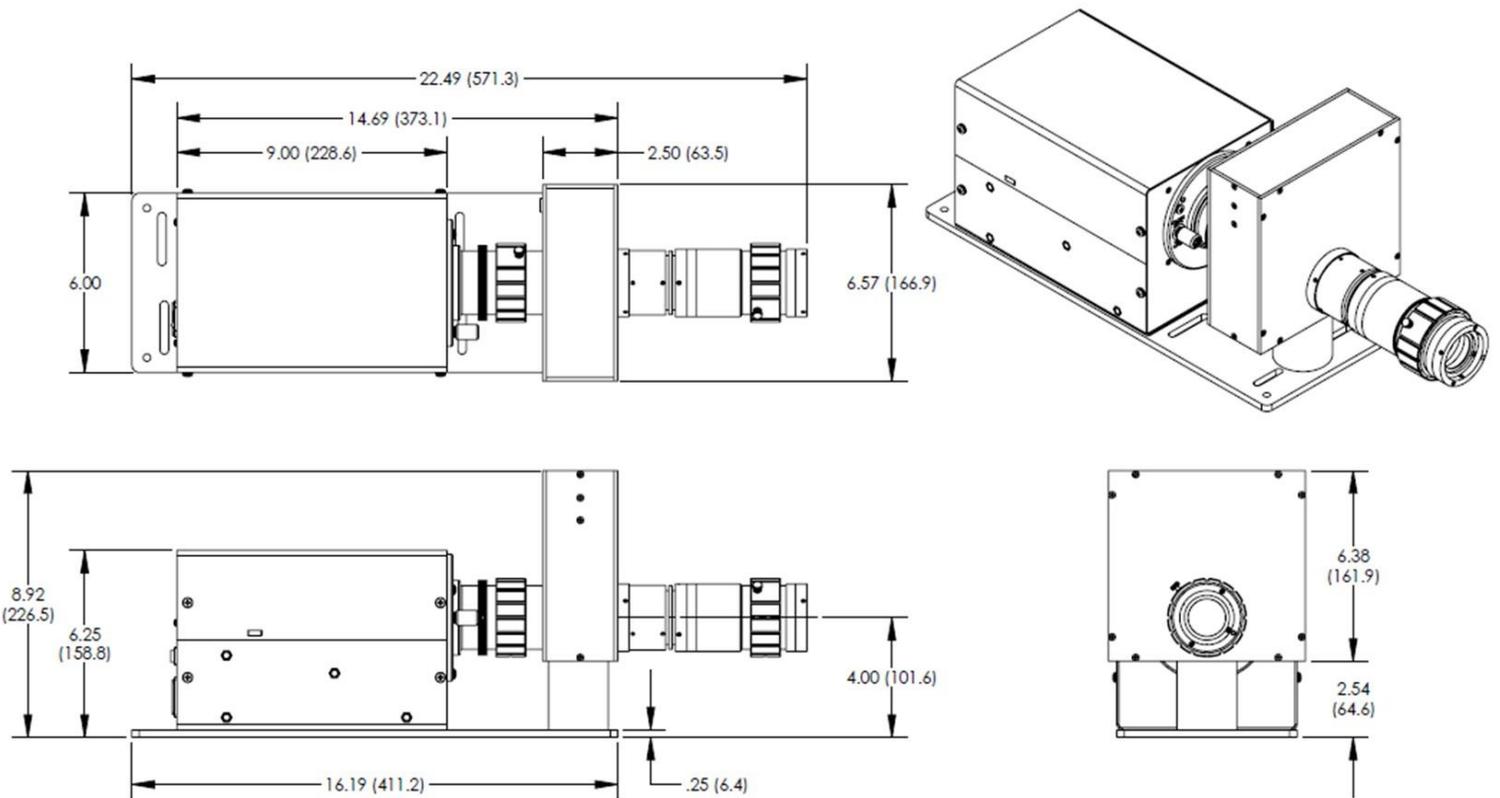


Figure 2: Dimensional Diagram

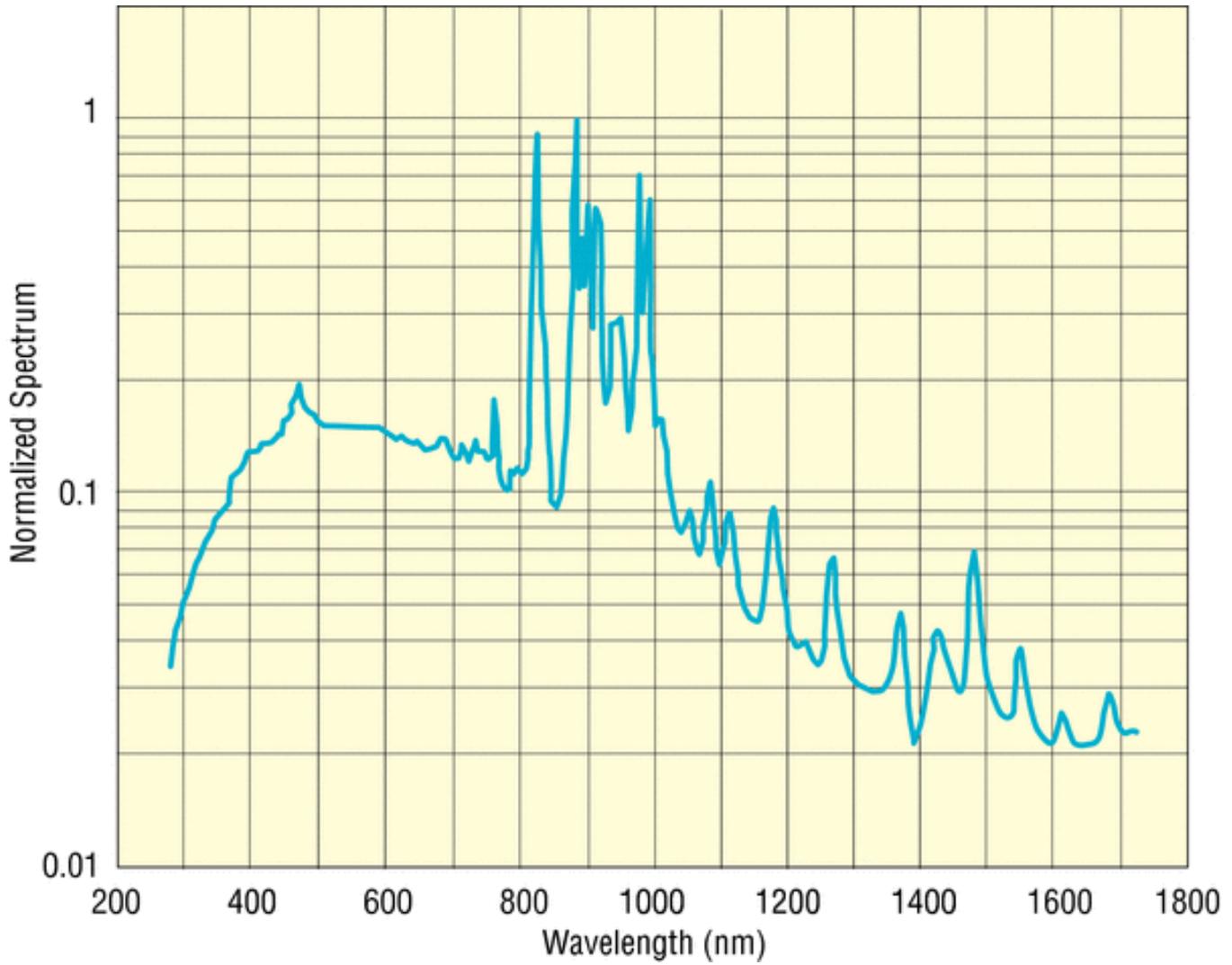


Figure 3: APEX2-XELamp Spectra

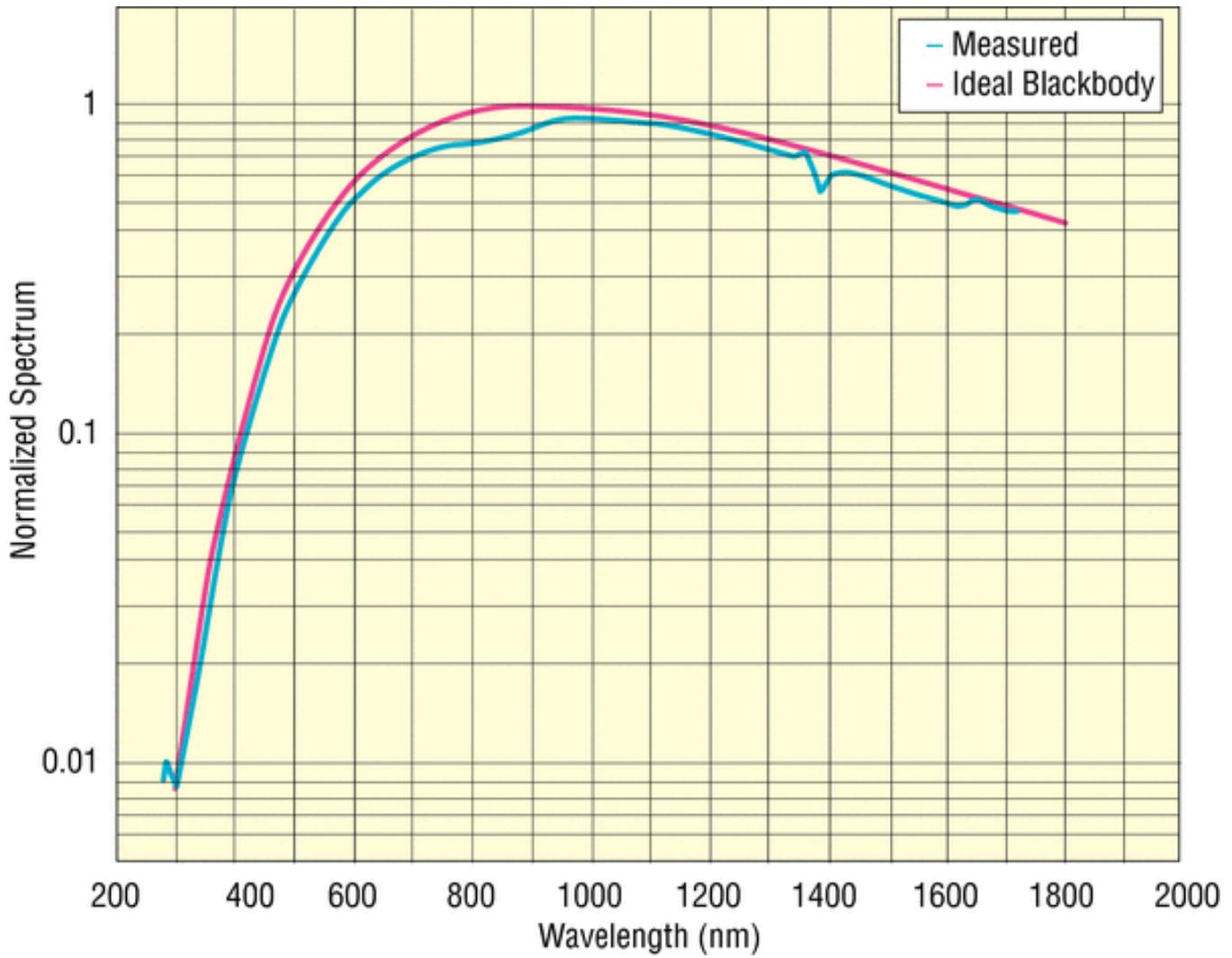


Figure 4: APEX2-QTH Lamp Spectra

3 SYSTEM SETUP

3.1 ITEMS INCLUDED WITH SYSTEM

Oriel Instruments provides a pre-aligned light source and filter wheel installed on a mounting plate. In addition, the system includes:

- 100 W Xe or QTH Lamp
- Front lens assembly
- Integrated arc lamp or QTH source
- USB stick preloaded with Universal Filter Wheel Utility Software, this User's Manual, and Quick Start Guide
- Filter wrench
- 1/16 hex wrench
- US AC power cord
- EU AC power cord
- 74010-1020 Filter wheel cable assembly
- 3 pcs 6-32 x 3/16 socket set cone point set screws
- 2 pcs 0.8 A, 250 V 5 x 20 m fuse
- 4 pcs mounting feet

3.2 UNPACKING

Remove all items from the shipping containers and verify each item is accounted for. The system is carefully packaged to minimize the possibility of damage during shipment. Inspect the shipping boxes for external signs of damage or mishandling. Inspect the contents for damage.

If any item is missing or damaged, immediately contact Oriel Instruments or the Newport representative from whom the system was purchased.

It is suggested to save the packaging material and shipping container, in case the equipment needs to be relocated at a future date.

	<p>WARNING</p> <p>Do not attempt to operate this equipment if there is evidence of shipping damage or there is suspicion that the equipment will not operate correctly. Damaged equipment may present hazards.</p>
---	---

3.3 CHOOSING A LOCATION

Choose an installation location where the power requirements can be met for the system. Be sure power is not applied to the system until the setup has been completed.

The environment should be that of a typical laboratory atmosphere, without excessive humidity and contaminants in the air. Do not allow the ventilation holes on the illuminator or its computer to be blocked. Air should be able to circulate freely around the system.

Carry the instrument by its base plate. When the system is placed in its final location, check to ensure that none of the pre-assembled items have become loose during transportation and handling.

3.4 ASSEMBLY

To ensure the front lens assembly and integrated light source components of the APEX2 unit remain undamaged during shipping, these two components are packaged as separate items within the protective foam the APEX2 is shipped in. Some minor assembly is required before the lamp can be installed.

1. The top piece of foam houses the miscellaneous accessories included with the APEX2 (AC power cords, etc.), front lens assembly, and Xe/QTH lamp. Remove the protective foam placed over the front lens assembly, remove the front lens assembly from its protective packaging, and set the front lens assembly aside.



Front Lens Assembly



Figure 5: How the APEX2 unit arrives in its packaging

2. Remove the top piece of foam to reveal the integrated arc lamp/QTH source. Remove the integrated light source from the protective foam. The two items as shown in Figure 6 are now ready for assembly.

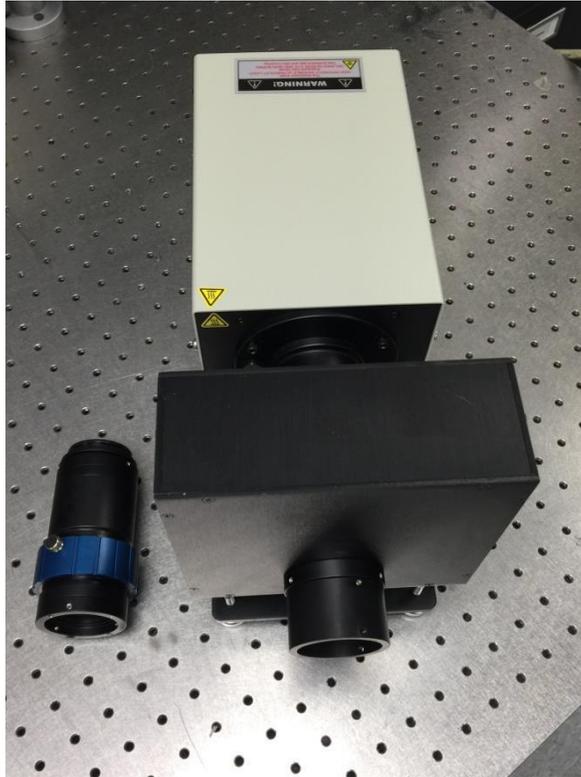


Figure 6: Front lens assembly and integrated light source

3. Insert the male flange of the front lens assembly into the female flange of filter wheel of the integrated light source unit as shown in Figure 7. Install with the set screw of the blue Collimating Lens Adjustment Ring oriented upright as shown in Figure 8. This will make the set screw easily accessible for future adjustment. When the front lens assembly is installed in an optimal orientation, secure the assembly in place with the three 6-32 x 3/16 socket set cone point set screws provided.

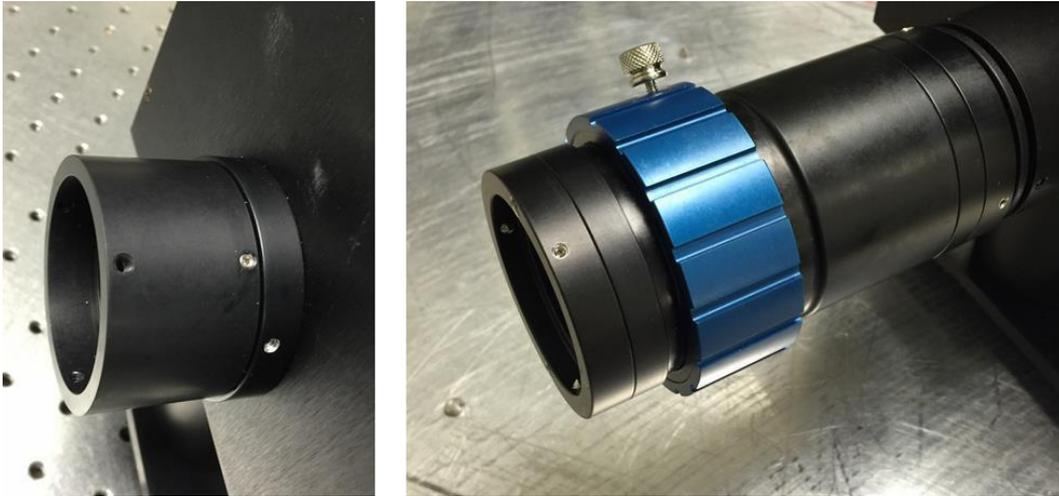


Figure 7: Female flange of the integrated light source unit (left) and installed front lens assembly (right)

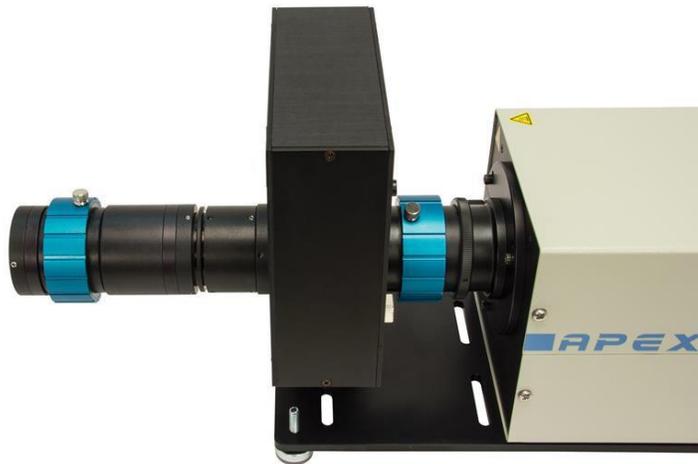


Figure 8: Installed front lens assembly with Collimating Lens Adjustment Ring set screw oriented vertically for easy accessibility

The APEX2 is now ready for lamp installation.

3.5 LAMP INSTALLATION

Before plugging in the power cord and powering up the APEX2, the lamp must be installed. The lamp is shipped in its own box to prevent damage during transportation. It is suggested to keep this box, as it can be utilized if moving or transporting the APEX2 in the future. Never move the APEX2 while the lamp is installed. Damage to the lamp or system may result.

	ELECTRICAL SHOCK HAZARD Do not open the APEX2 cover and attempt to work inside without first turning the instrument off and disconnecting the power cord from the AC mains. Failure to follow this warning can result in severe injury or death.
---	--

	WARNING When installing the lamp, it is <u>required</u> to: <ul style="list-style-type: none">• Wear eye protection.• Wear powder-free gloves.• Make sure the power is turned off.
---	--

1. Using a Phillips head screwdriver, remove the (4) screws holding the top cover of the housing and set aside. Refer to Figure 9.



Figure 9: Remove Top Cover Screws

2. Remove the top cover and set aside as shown in Figure 10.

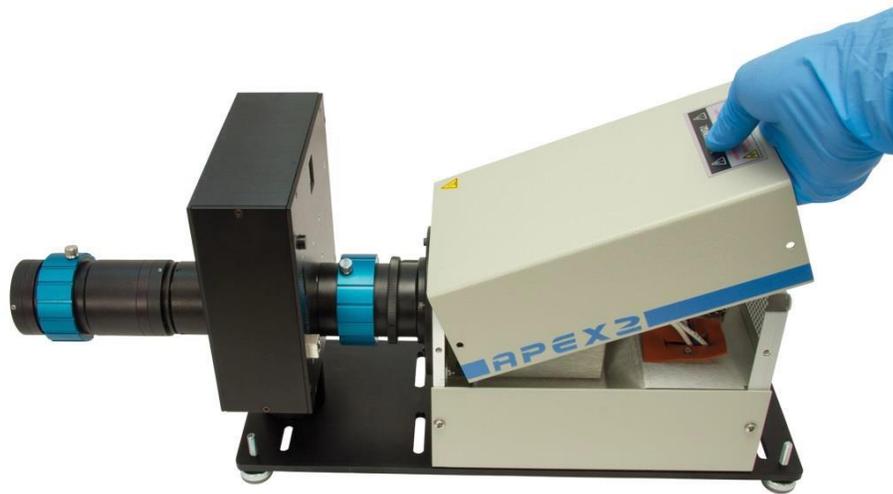


Figure 10: Remove Housing Cover

3. Loosen and remove the knurled thumb screw as shown in Figure 11. Remove the lamp cable holder plate and set aside with the thumb screw.



Figure 11: Remove Lamp Cable Holder

4. Press the lamp holder wires toward each other and then pull them toward the rear of the APEX2 unit. This is demonstrated in Figure 12.

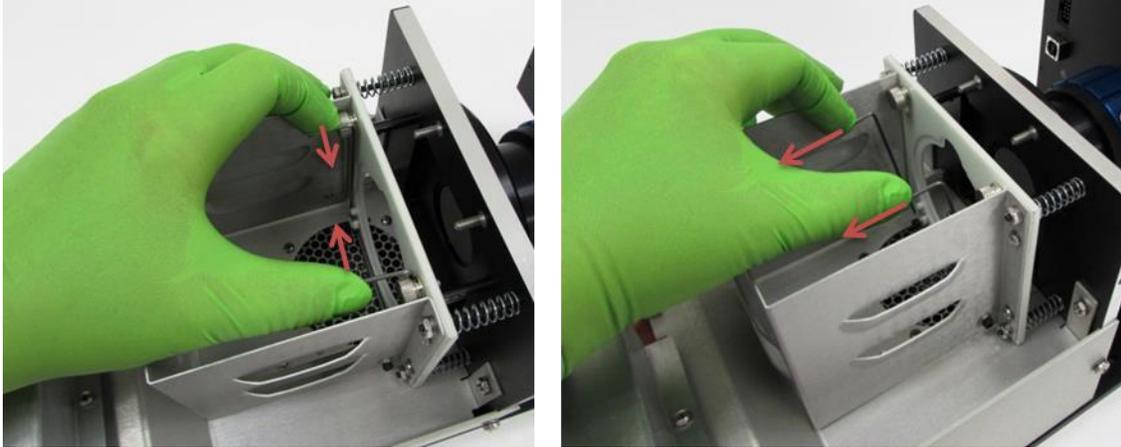


Figure 12: Move Lamp Holder Wires

5. The power supply cable shown in Figure 13 is secured for transportation. Unfasten it and let it hang to the side for now.

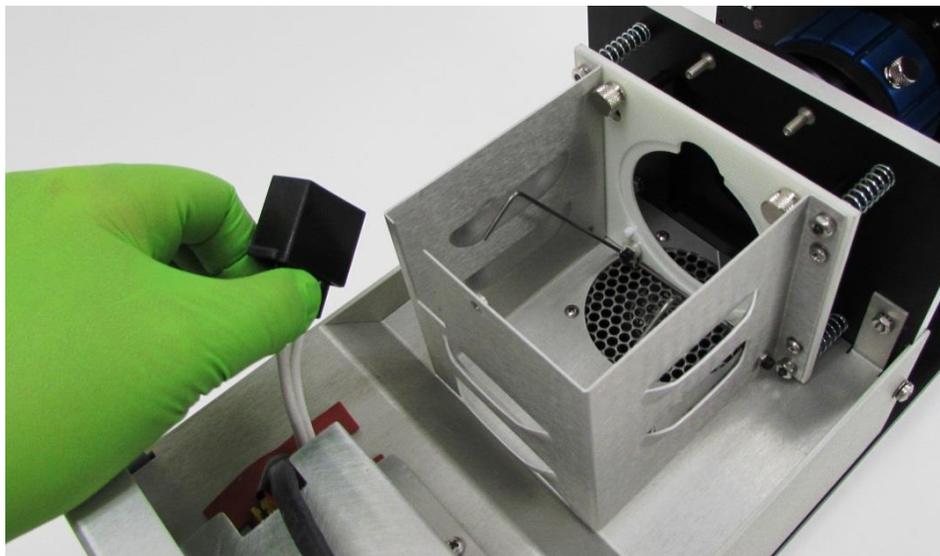


Figure 13: Unfasten Power Supply Cable

6. Insert the lamp as shown in Figure 14. The outer edge of the elliptical reflector must fit into the groove shown in Figure 15.



CAUTION

Always wear powder-free gloves or finger cots when handling the lamp. Handle the lamp from the outside of the reflector. Making contact with the lamp or reflector may cause permanent damage.



Figure 14: Insert Lamp

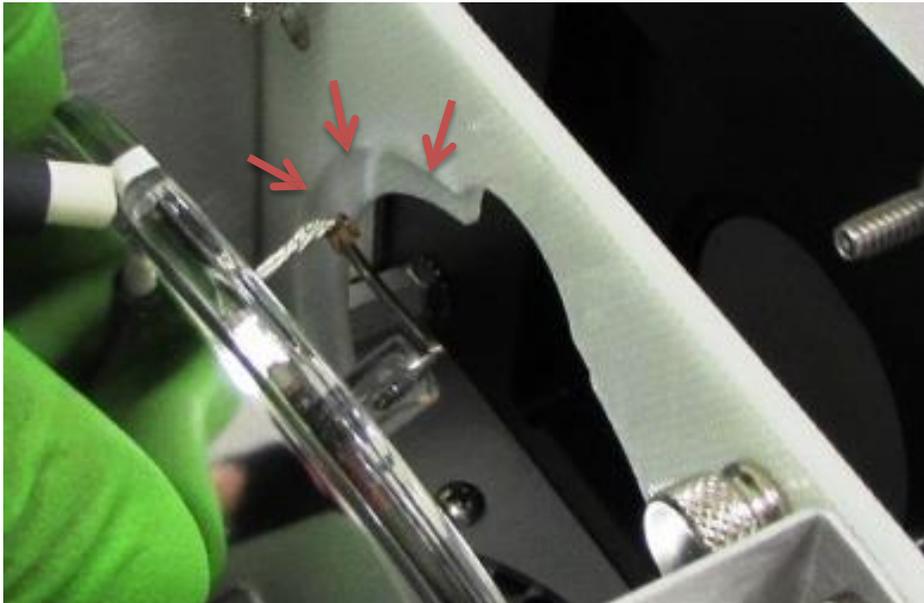


Figure 15: Lamp Ellipse Positioning Groove

7. Secure the lamp in place by returning the lamp holder wires to their original location per Figure 16.

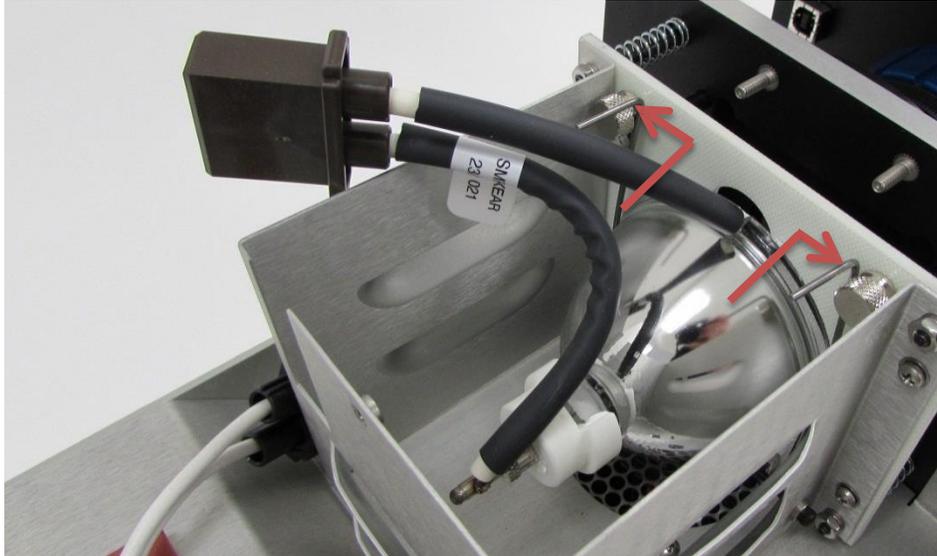


Figure 16: Secure Lamp

8. Connect the lamp wiring to the power supply cable. The connectors are designed so that they cannot be connected incorrectly, so there is no need to be concerned with lamp polarity. See Figure 17.



Figure 17: Connect Lamp Wiring

9. Re-attach the lamp cable holder which was set aside earlier by tightening the thumb screw as shown in Figure 18. The completed lamp installation is shown in Figure 19.

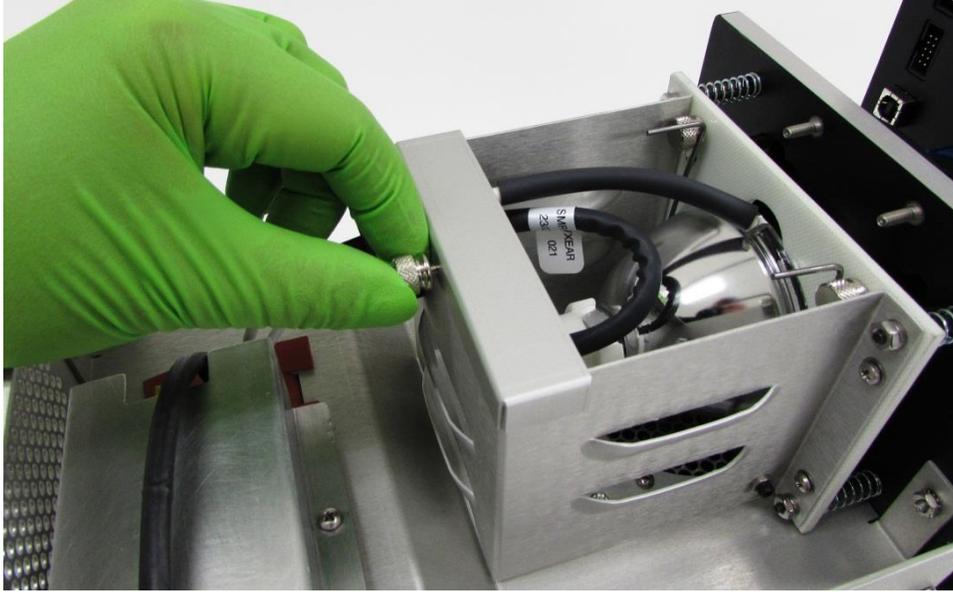


Figure 18: Re-attach Lamp Cable Holder

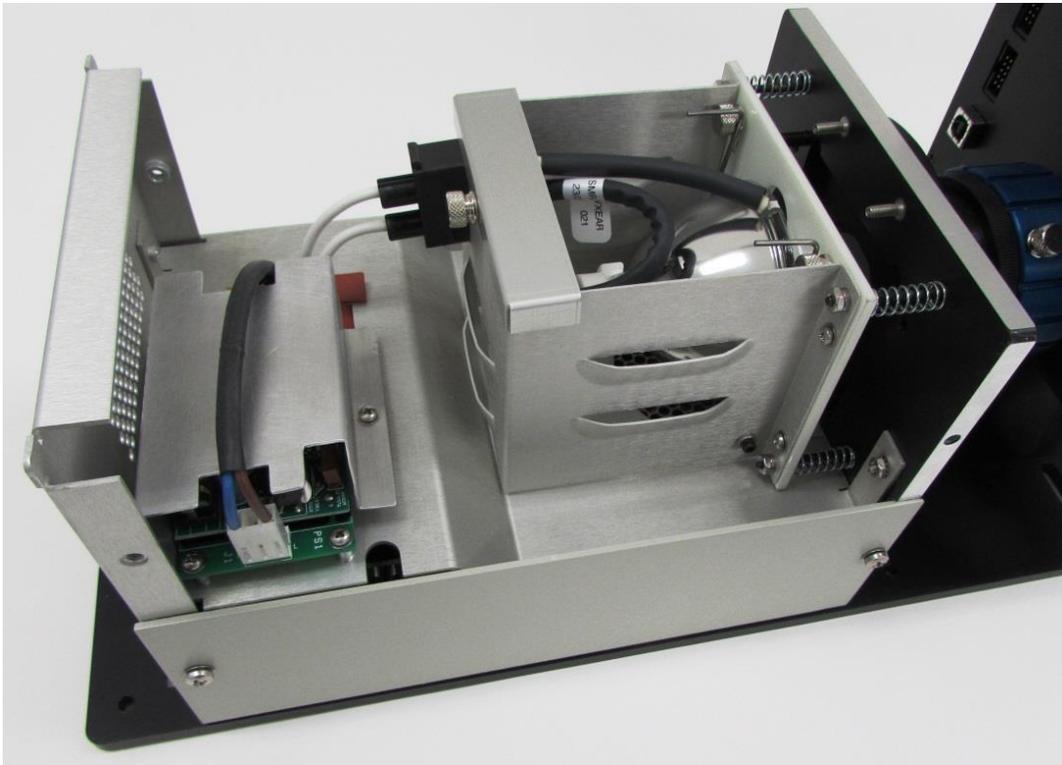


Figure 19: Completed Lamp Installation

10. Re-install the cover, ensuring that the top cover's safety interlock lines up with its slot in the bottom portion of the housing as shown in Figure 20.



Figure 20: Safety Interlock

11. Using a Phillips head screwdriver, re-install the (4) screws to secure the top cover of the housing. Do not leave out any screws.

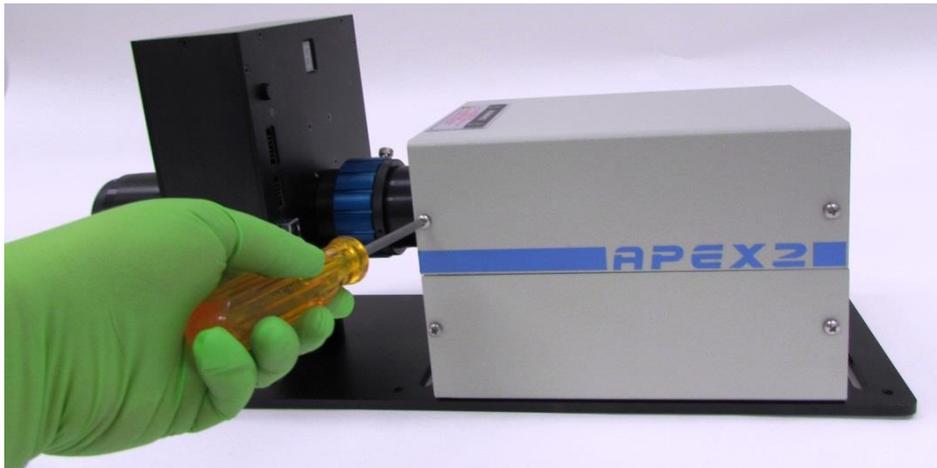


Figure 21: Cover Re-installed

3.6 ELECTRICAL CONNECTION

Before powering up the system for the first time, it is suggested to have a qualified electrician verify the wall socket to be used with the APEX2 meets the requirements for operation as noted.

The line voltage requirements for the APEX2 are as follows:

100 to 120 VAC, 50-60 Hz
200 to 240 VAC, 50-60 Hz

	<p>ELECTRICAL SHOCK HAZARDS</p> <p>To avoid electric shock, connect the instrument to a properly earth-grounded, 3-prong receptacle only. Failure to observe this precaution can result in severe injury or death.</p> <p>Never attempt to open the lamp power supply or monochromator power adapter. These items do not contain any user serviceable parts. Failure to follow this warning can result in severe injury or death.</p>
---	--

The APEX2 illumination systems conform to CE standards for both safety and EMC. During normal use, this equipment will not pose any electrical hazards to the user. Read all warnings before installing or operating this system. Never open the APEX2 cover and attempt to work inside without first turning the instrument off and disconnecting the power cord from the AC mains.

	<p>CAUTION</p> <p>Do not turn on the APEX2 until all connections have been made to the monochromator, computer, hand controller and/or external 5 VDC source. Refer to Section 4 for information on various methods and required connections to control the APEX2.</p>
---	---

If there are any questions or concerns, contact Oriel Instruments or the regional sales representative for Newport.

	<p>WARNING</p> <p>When the power switch is turned on, the lamp will begin emitting light immediately. Do not turn on the APEX2 until the output flange is directed in such a way that people, animals and equipment will not be harmed by the light.</p>
---	---

A dedicated power line or line isolation may be required in certain locations, as the electronics contained in the instrument are sensitive to static electricity and radiated electromagnetic fields. Operation of the system near intense pulsed sources (lasers, Xenon flash lamps, etc.) may compromise performance.

Before making any electrical connections, verify the power switches are in the OFF position for the APEX2 and monochromator, if being used.

Connect the power cord to the back of the APEX2 before to plugging it into an AC electrical outlet.



Figure 22: APEX2 Power Connection

NOTE: Always allow the lamp to warm up for at least 30 minutes prior to beginning work.

3.7 LAMP INTENSITY ADJUSTMENT (APEX2-QTH ONLY)

The model APEX2-QTH allows for lamp intensity to be controlled by an external DC voltage source. External control is enabled by moving the toggle switch on the back panel to the **UP** position. When the toggle switch is in the **DOWN** position, the lamp will operate at its full power rating of 100 watts.

The BNC female connector on the back of the APEX2-QTH is used to connect a DC voltage source. The input voltage range is 0 to 5 VDC. 0 VDC corresponds to the lamp being off; 5 VDC allows the lamp to run at its maximum power of 100 watts.

NOTE:

In the event that the lamp is not operating and an external voltage source is not being applied, check the position of the toggle switch and ensure it is in the DOWN position.

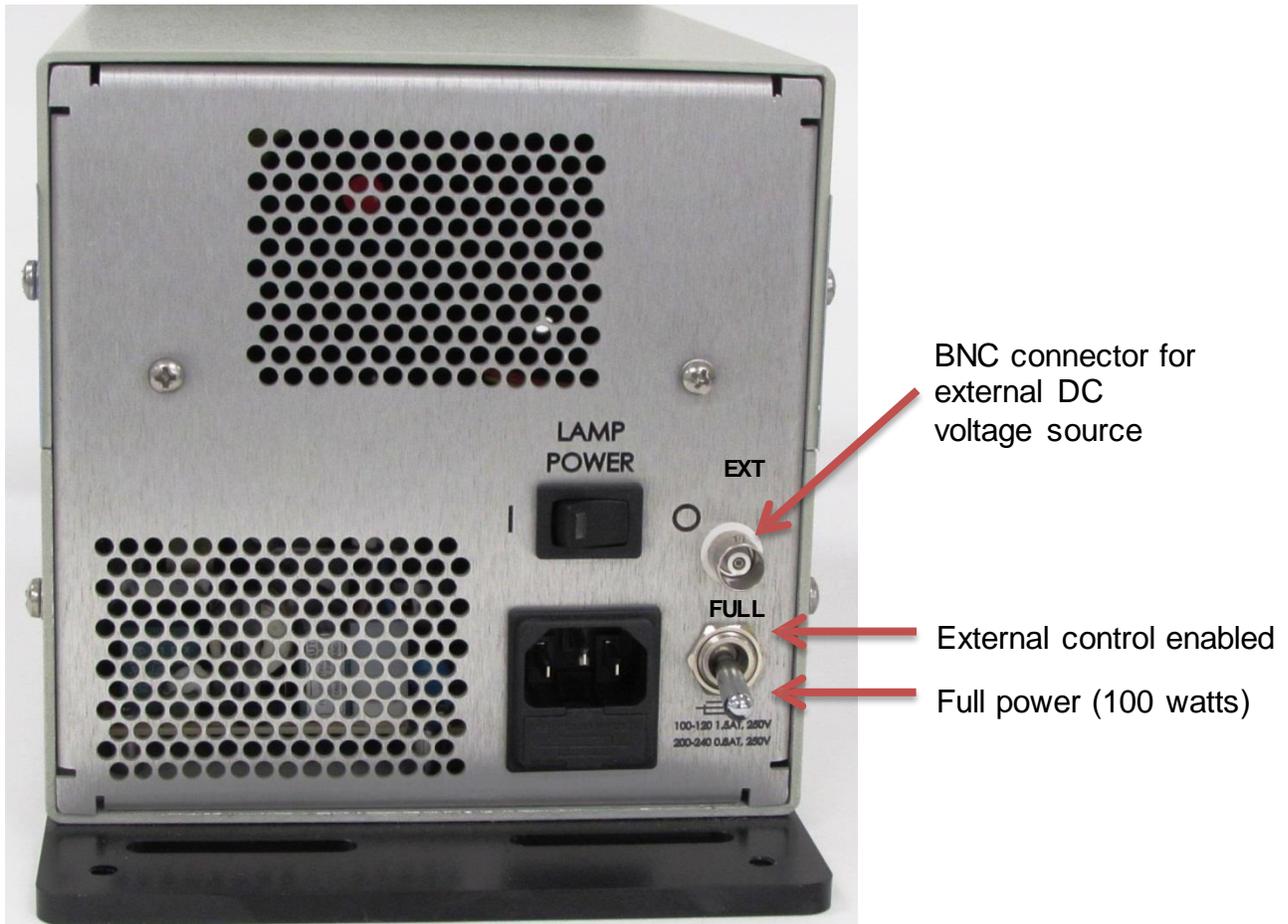


Figure 23: Lamp Intensity Adjustment Settings

4 CONTROLLING THE APEX2 FILTER WHEEL

The APEX2 integrates a light source, filter wheel and power supply into a complete package. Power is supplied to both the lamp and the filter wheel. A large, easy to read digital display indicates which filter is active, i.e. which filter is in the light path. The digital display is positioned on the rear panel of the filter wheel assembly, so that it may be observed from a safe and convenient viewing angle not within the path of the output light.

The filters may be selected using the following:

- Manual pushbutton control
- Direct APEX2 computer control
- Oriel Monochromator Utility program
- Monochromator command set
- Monochromator Hand Controller
- Oriel TracQ Basic radiometry software

For those who wish to create their own programs, low level commands may be sent to the system via its USB connection to a PC. Filter installation is covered in Section 6.

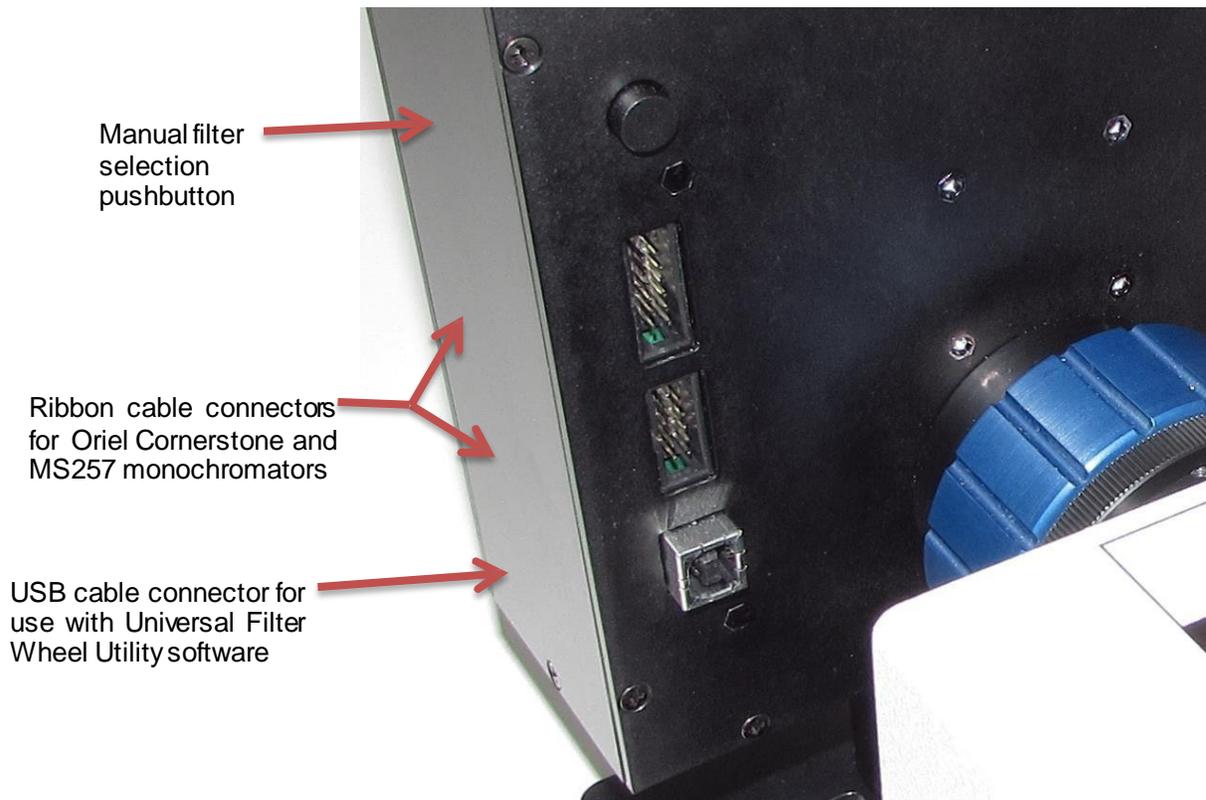


Figure 24: Filter Wheel Connectors

4.1 MANUAL CONTROL

Pressing the pushbutton located on the back of the filter wheel, as shown in Figure 25, advances the filter by one position.

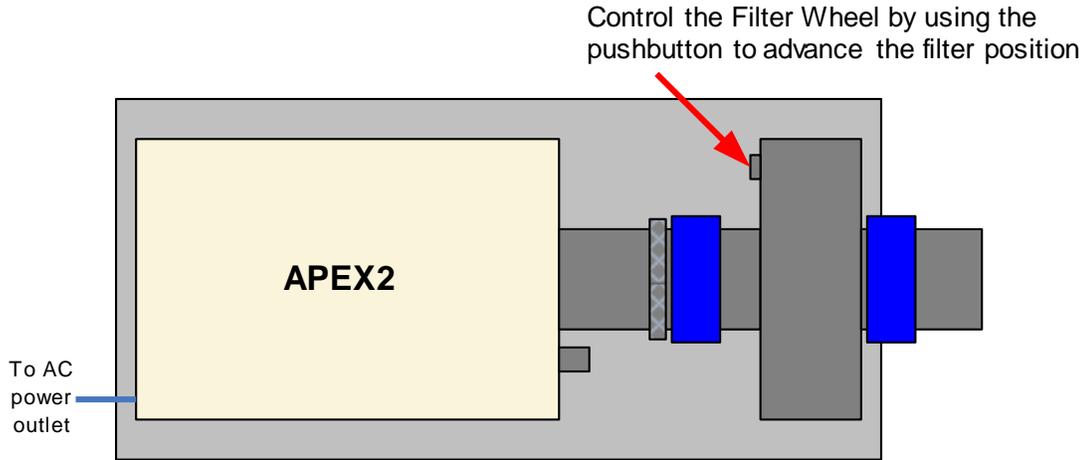


Figure 25: Manual Control Configuration

4.2 DIRECT APEX2 COMPUTER CONTROL

The APEX2 filter wheel comes with its own utility program and command set. To utilize these control options, connect the filter wheel directly to a computer using a USB cable. Refer to Figure 26 for wiring connection information. Section 7 discusses the utility program and command set in detail.

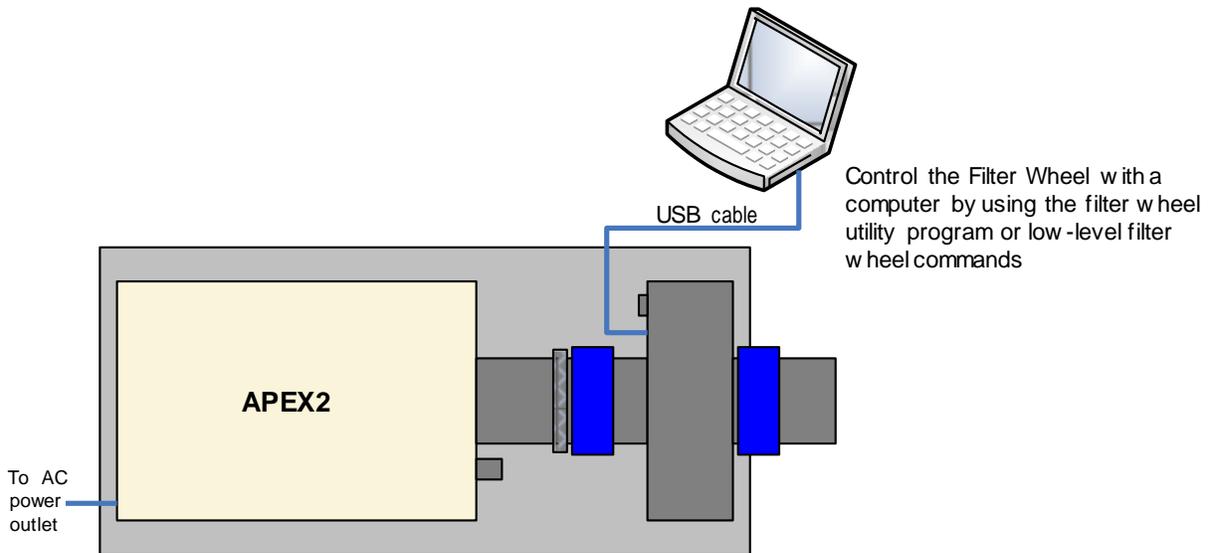


Figure 26: Direct Computer Control Configuration

4.3 MONOCHROMATOR UTILITY SOFTWARE CONTROL

Oriel's Cornerstone and MS257 series monochromators include a utility program used to query and command the instruments. When the filter wheel is connected to the monochromator, the monochromator's utility program may be used to control both instruments. Refer to the monochromator user's manual for more information.

Ensure the power switches switch for the APEX2 and monochromator are in the off position (marked as O). Then connect the power adapter and all other cables to the monochromator, as shown in Figure 29.

The monochromator is connected to the APEX2 filter wheel using a ribbon cable provided with the APEX2 monochromator adapter kits. Adapter kits are listed in Section 8.1. Connect the ribbon cable to the filter wheel using whichever connector matches the pin configuration of the cable.

Connect the other end of the power cord to the AC mains only after the lamp has been installed and all other connections made. Once the APEX2 and monochromator are powered on, the monochromator allows the user to select which filter is placed in the optical path.

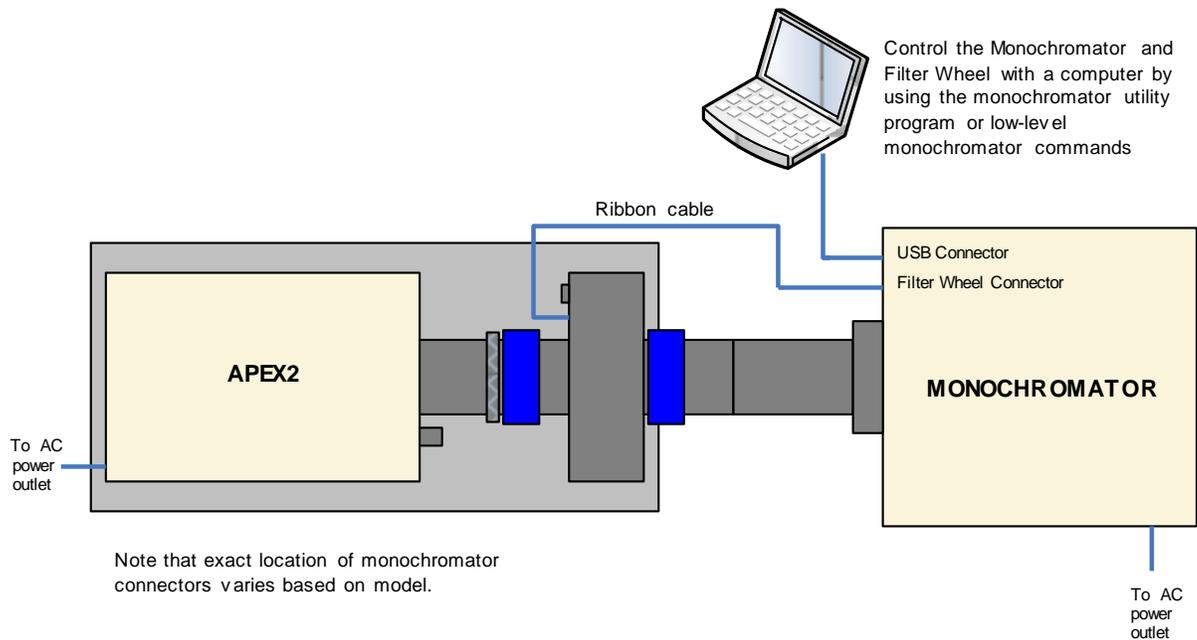


Figure 27: Monochromator Control Configuration

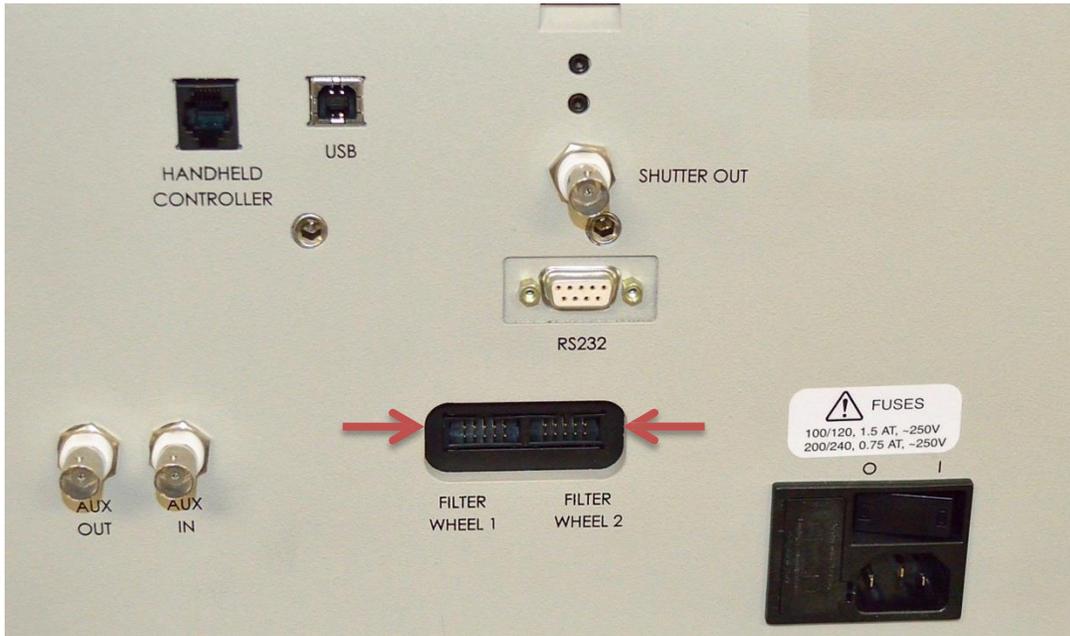


Figure 28: MS257 Monochromator Connections

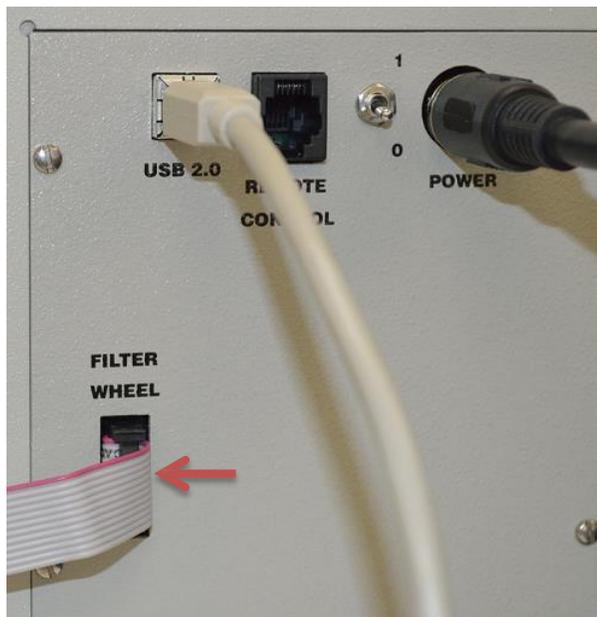


Figure 29: Cornerstone USB Monochromator Connections

4.4 MONOCHROMATOR COMMAND SET

For those who want to incorporate monochromator as well as APEX2 filter wheel control into their own custom software, both instruments may be controlled by the command set used by the monochromator.

Set up the monochromator and APEX2 as noted in Section 4.3 and Figure 27. Refer to the monochromator's user manual for information on the command set.

4.5 MONOCHROMATOR HAND CONTROLLER

For those who do not wish to use a computer to control the Cornerstone Series and MS257 Series monochromators, Oriel offers hand controllers which plug directly into the monochromators. When the APEX2 is connected to the monochromator via a filter wheel ribbon cable, the hand controller can be used to select a filter. Refer to the monochromator's user manual for more information.

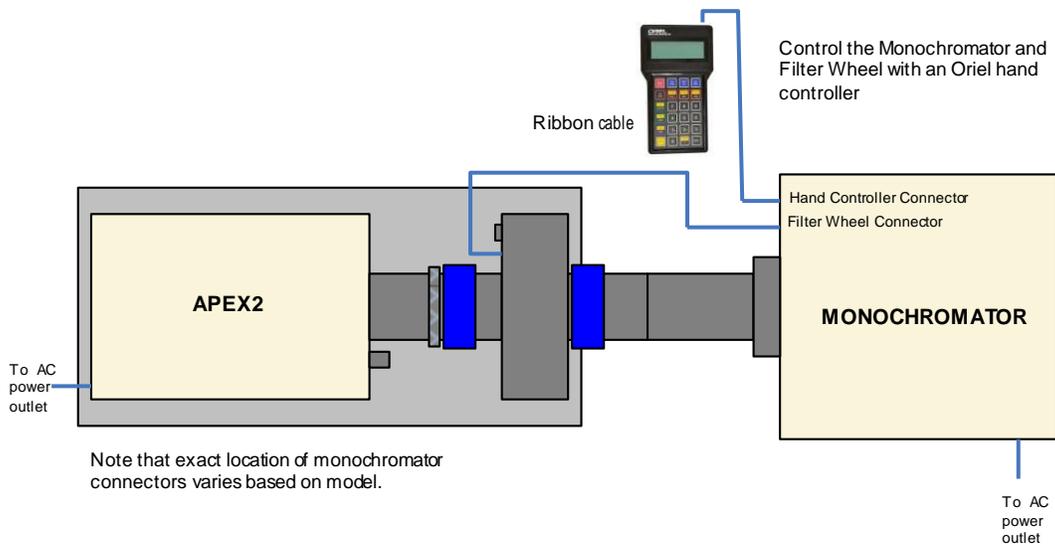


Figure 30: APEX2 Configured with Hand Controller

4.6 TRACQ BASIC CONTROL

In order to ensure that the light exiting the monochromator is truly monochromatic, it is necessary to incorporate optical filters into the design. The physics of diffraction gratings is such that higher order wavelengths need to be blocked. The filters installed into the APEX2 need to be chosen to block unwanted wavelengths.

By setting the filter parameters in the TracQ Basic software, the monochromator automatically selects the appropriate filter for any desired wavelength. Refer to the TracQ Basic manual for more information.

Figure 31 shows the filter setup screen for a Cornerstone monochromator. The MS257 setup screen is similar, but uses five filters and up to two filter wheels simultaneously.

Filter Selection

Select Filter

Present System Status

Filter In Use	Filter Label
2	Using Table
Current Wavelength	Wavelength Units
555.00	nanometer

Automatic Filter Change-Over Table

From filter	to filter	at wavelength
1	2	355.00
2	3	625.00
3	4	1100.00
4	5	1580.00
5	6	Inf

Figure 31: Filter Selection Table in TracQ Basic

NOTE:

The parameters entered in Figure 31 are an example. The actual filter changeover parameters depend on the monochromator configuration and filters being used.

5 APEX2 FEATURES

The APEX2 provides a number of features into a convenient, compact illumination system. All features are integrated into the unit, to keep the system easy to transport, set up and operate.

5.1 FILTER WHEEL OVERVIEW

The motorized filter wheel holds up to six (6) filters. The MS257 monochromator was designed to work with filter wheels containing up to five (5) filters. When the MS257 is used with the APEX2, the first five filters will be selectable. When used alone or with other monochromators, all six (6) filter positions may be utilized. The active filter position change at a rate of 1.7 seconds per filter. When applying power to the APEX2, Refer to Section 4 regarding the various methods by which the filter wheel may be controlled and Section 6 for filter installation.

5.2 SHUTTER

The integrated shutter is a very convenient feature which may be used to close off the light path while the lamp is warming up, or a change in setup is being performed. Frequently restarting the lamp may shorten the lamp's life and resets the warm-up time. If the lamp is not needed for a short amount of time, the shutter can simply be closed.

Turning the knob clockwise opens the shutter. Rotating the knob counterclockwise closes the shutter.

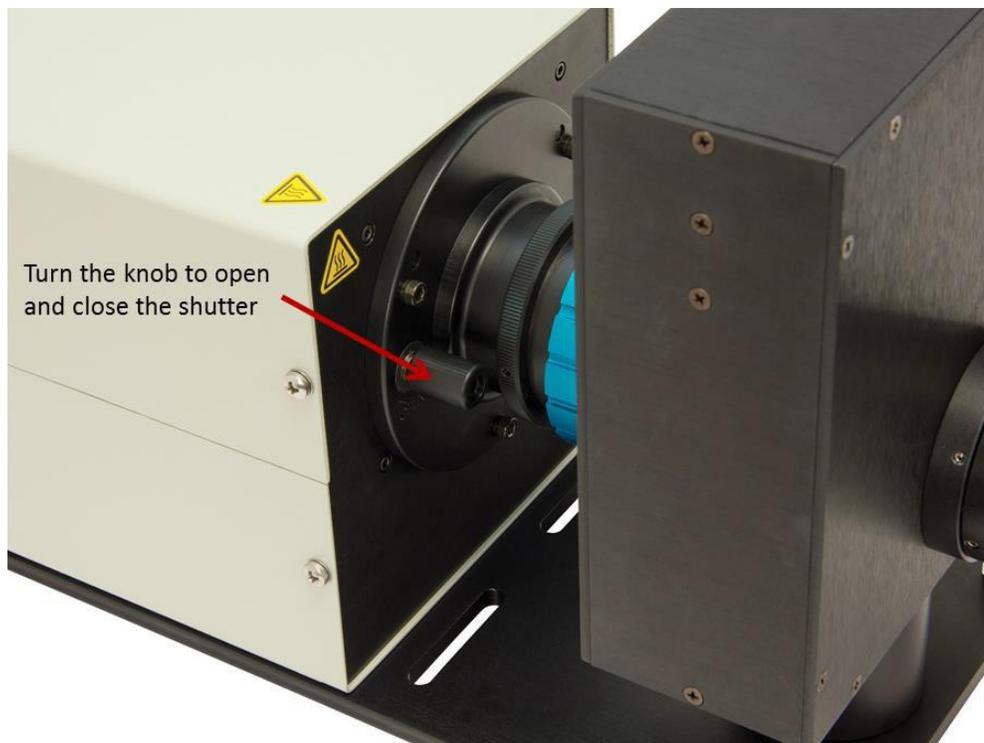


Figure 32: Shutter Knob

5.3 ADJUSTABLE COLLIMATING LENS

The collimating lens may be adjusted to optimize light coupling. Loosen the silver thumb screw shown in Figure 33. Turn the blue knob near the output flange to change the lens position. Be sure to tighten the locking screw when finished.

Note: under normal circumstances, this is the only lens which should be adjusted. Do not make any changes to the relay lens position, as its position was optimized before shipment.



Figure 33: Collimating Lens Adjustment

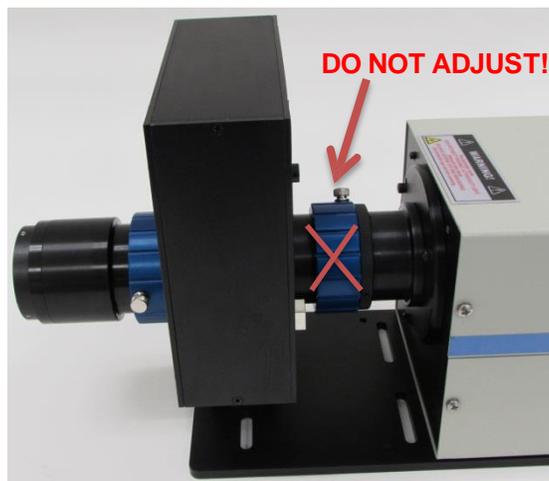


Figure 34: Do Not Adjust Relay Lens

5.4 MATING FLANGE OPTIONS

The APEX2 is designed with a standard Oriel 1.5-inch series mating flange, allowing it to be coupled to a variety of instruments and accessories. Items are connected to the mating flange by tightening three set screws, shown in Figure 35.



Figure 35: Oriel 1.5-Inch Series Flange

6 FILTER INSTALLATION

	<p>WARNING</p> <p>When the light source is in use, the filter wheel and all other items in the optical path may become hot. Ensure items are cool before handling them.</p> <p>Ensure the shutter is closed and the lamp is off whenever working in the filter wheel compartment.</p>
---	--

Remove the front plate of the filter wheel assembly by unscrewing eight (8) flat head screws using a Phillips screwdriver. Refer to Figure 36.

Set all items aside, ensuring that no foreign matter can contaminate or damage the optics. Do not attempt to disconnect the output optics and flange from the front plate.



Figure 36: Removing the Front Panel of the Filter Wheel

Each filter position is labeled on the wheel. Remove the filter retaining ring using the filter wrench included with the system. Insert the filter and re-install the retaining ring. Always wear gloves or finger cots when handling optics. See Figure 37 and Figure 38.

For all filter positions which are not being utilized, ensure the retaining rings are snug so that they do not become loose and obstruct the filter wheel. Only then should the filter cover be re-installed.

Note: Always keep track of which filters are installed at each slot in the filter wheel. It is suggested to write down this information and update it whenever the configuration is changed. This is particularly helpful when multiple users will be operating the APEX2 system.



Figure 37: Filter Wrench Removing Retaining Ring

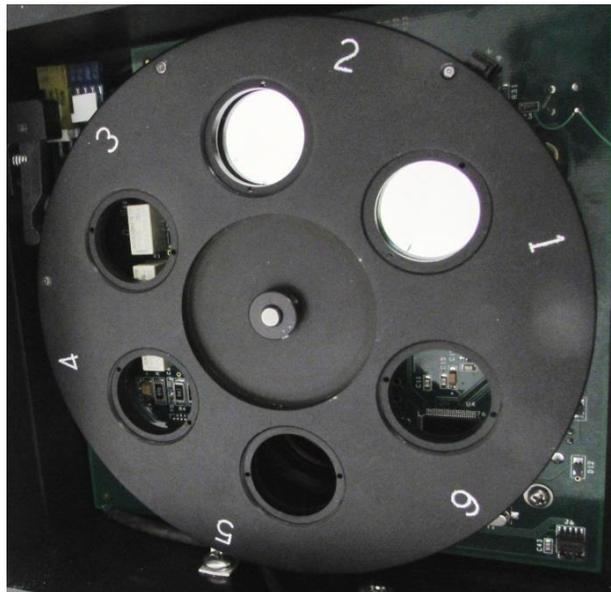


Figure 38: Installed Filters

7 UNIVERSAL FILTER WHEEL UTILITY SOFTWARE

The following procedure acquaints a first time user with the installation and use of the APEX2 utility software. Do not be concerned if your display does not exactly agree with the figures shown. The focus of this manual is to help the first time user learn how to install and use the software. It is highly recommended that the first time user read these instructions before attempting to use or install the software. To install the software, the APEX2 utility software, a compatible computer with a USB port, and an APEX2 illuminator are required.

Installation of APEX2 utility program will create:

- a folder called C:\Program Files\Newport - Oriel Instruments\Universal USB Filter Wheel (an alternative directory may be chosen during the installation process)
- an icon called "USB Filter Wheel Utility" in the start menu
- an icon for NI Max on the desktop

7.1 SYSTEM REQUIREMENTS

Operating System	Microsoft Windows XP, Service Pack 3 Microsoft Windows 7, 32-bit and 64-bit Requires Microsoft .net Framework 4.0
Processor Speed	2 GHz minimum
Random Access Memory (RAM)	1 GB minimum
Peripheral Hardware	USB 2.0 port
Hard Drive Space	800 MB minimum

7.2 SOFTWARE INSTALLATION PROCEDURE

1. Turn on the computer following the manufacturer's instructions. Before installing the software, close all applications. Disconnect all instruments from the computer before installing this software.
2. Insert the USB stick included with the APEX2 system into an unused USB port on the PC. The "AutoPlay" dialog box may open. Select "Open folder to view files." If the "AutoPlay" does not start installing the software automatically, navigate to the USB stick's directory in Windows Explorer.
3. Double click the Setup.exe application, shown in Figure 39, application to begin the installation.

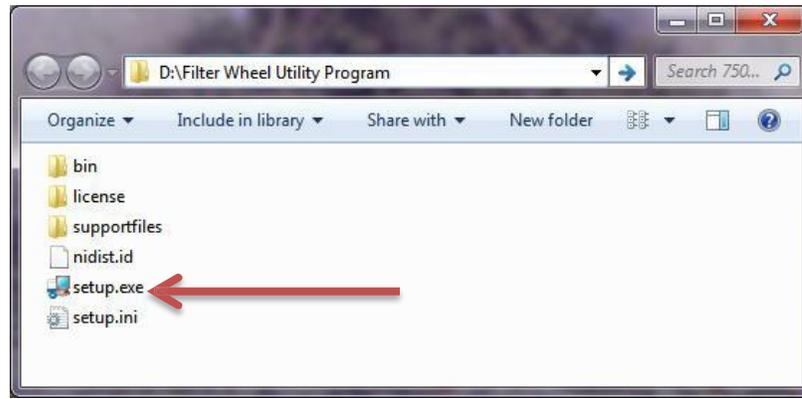


Figure 39: Setup Application

4. When the "Destination Directory" screen appears, the option is available to install the software into the default location or a different location. To install software into a different location, click the "Browse" button and select another directory. There are separate directories for APEX2 and National Instruments Products, as shown in Figure 40. Click "Next" to continue. During the installation, it is normal for other dialog boxes to flash as software is added.

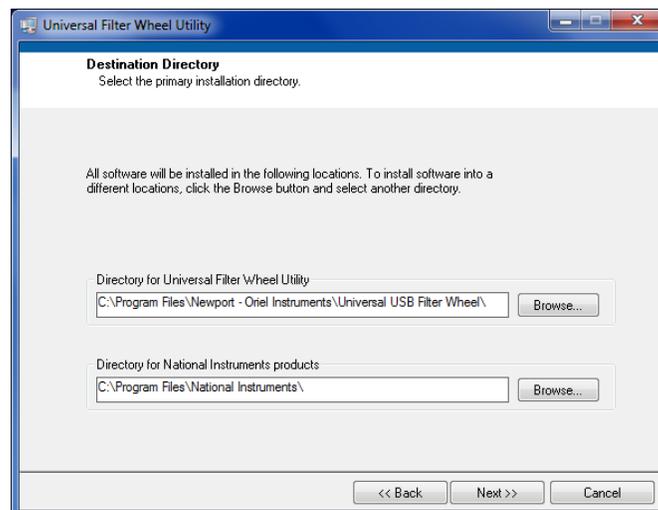


Figure 40: Destination Directory Selection

5. Choose to accept the license agreements and click Next.

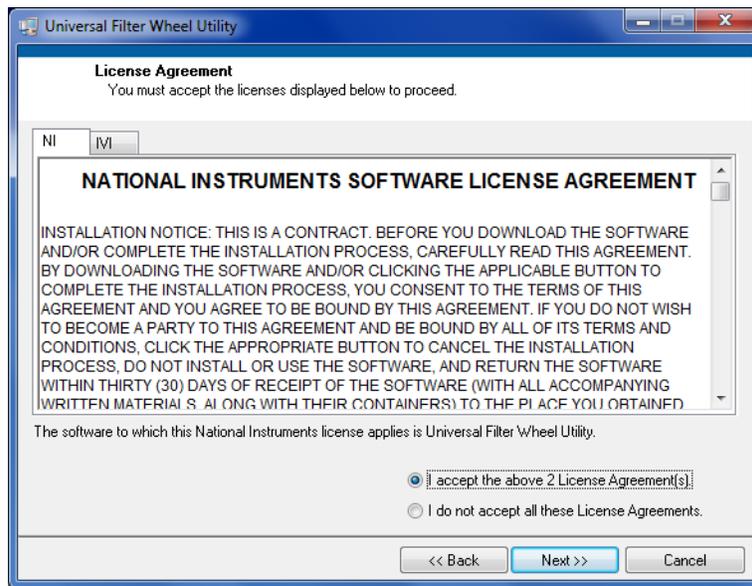


Figure 41: Accept License Agreements

6. Click "Next" to start installation. Otherwise click "Back" to change the installation settings.

Note: Clicking "Save File" will save a file in Rich Text File format (*.rtf). The .rtf file will list the installation summary text. This file will be named "save as install.rtf".

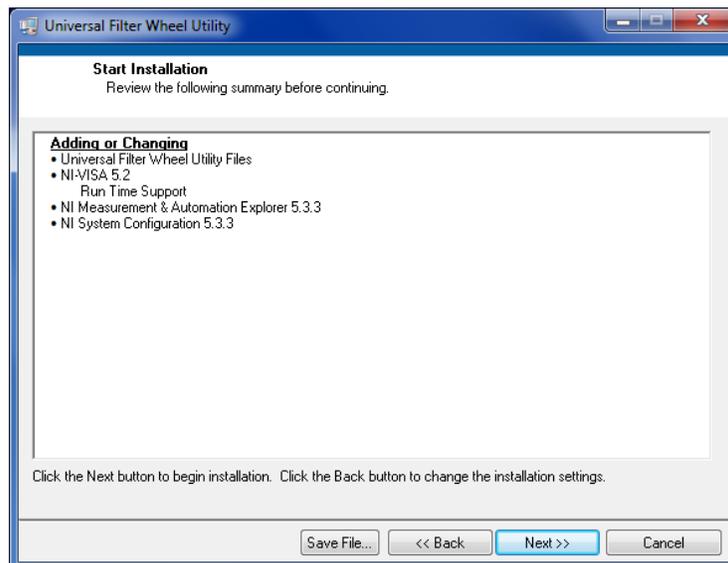


Figure 42: Start Installation

7. During installation, a status indicator will show the overall progress. Selecting "Cancel" brings up a dialog box to verify cancellation.

- When the installation is complete, click "Finish". Refer to Figure 43.

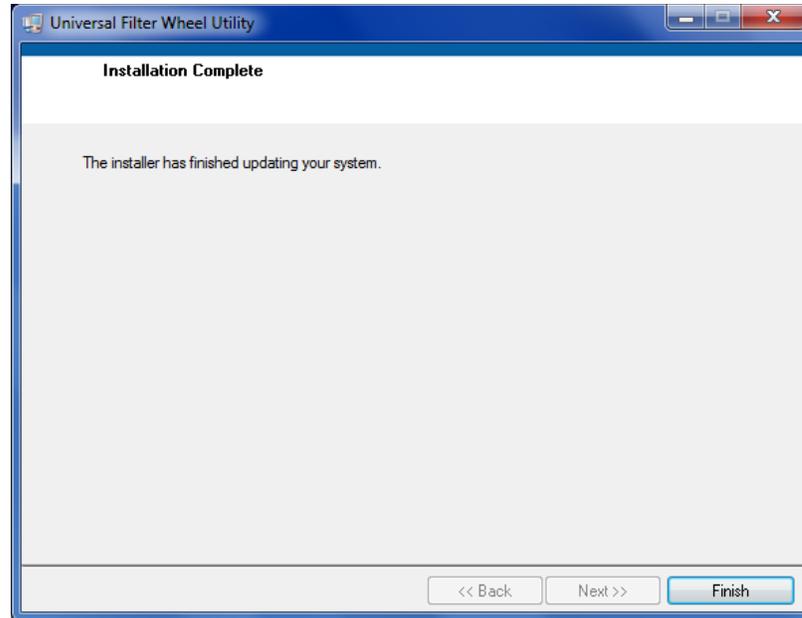


Figure 43: Finish Installation

- A dialog box will appear to restart the computer, as shown in Figure 44. This software can be used only after the computer is restarted.

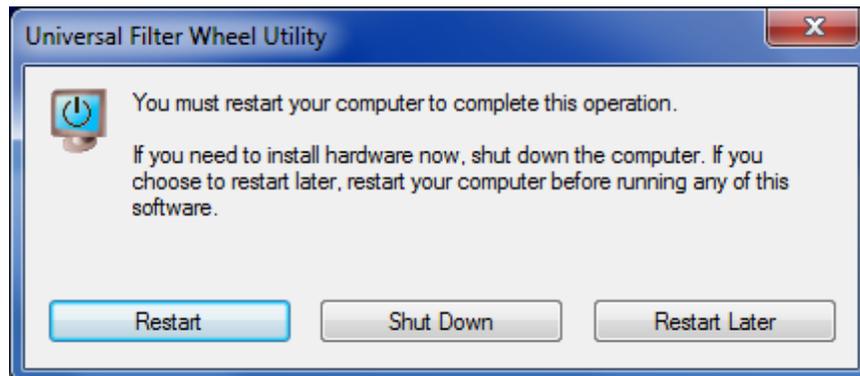


Figure 44: Restart Computer

- After the computer restarts, the Start menu shall contain the Filter Wheel Utility software and the National Instruments "NI Max" icon shall appear on the desktop, as shown in Figure 45.

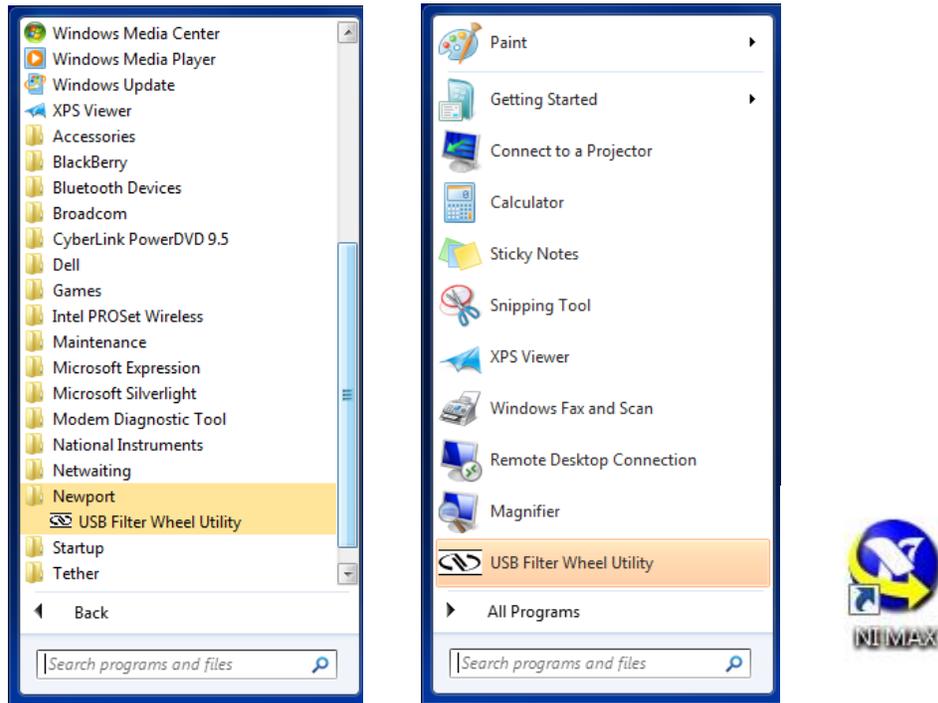


Figure 45: Start Menu and Desktop Changes

12. Right click the USB stick icon in the System Tray or Notification Area. When the pop-up menu shown in Figure 46 appears, click “Eject Mass Storage.” Wait for the “Safe To Remove Hardware” notification to appear as shown before removing the USB stick from the computer.

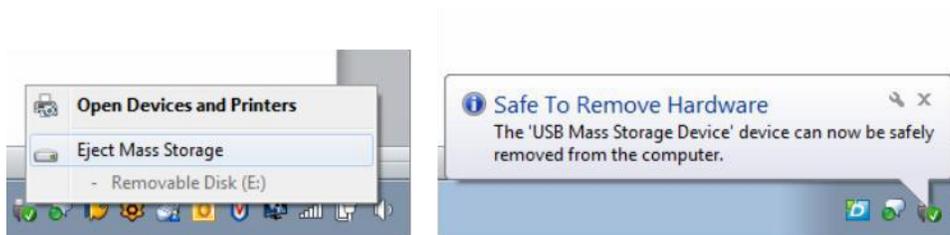


Figure 46: Safely removing the USB stick

7.3 USING THE FILTER WHEEL UTILITY SOFTWARE

1. After the Filter Wheel Utility software is installed and the computer restarted, connect the APEX2 to the computer via a USB cable and turn on the illuminator.
2. Run the utility program.
3. Select the instrument, and click the “Connect” button.

NOTE: If the software is unable to connect, check in the device manager to ensure that the USB driver has been installed correctly and the instrument is listed.



Figure 47: Connect to Filter Wheel

4. To verify communication, click “Device ID?”

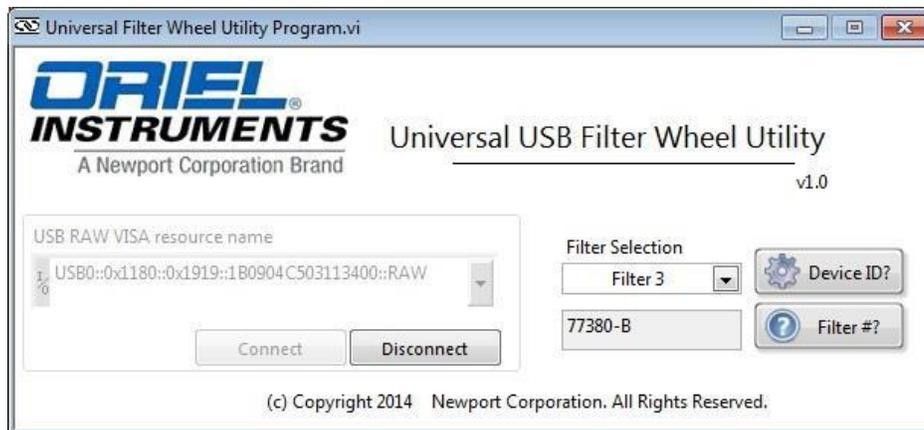


Figure 48: Device ID Query

5. To select a filter, click the pulldown menu and choose the desired filter position.

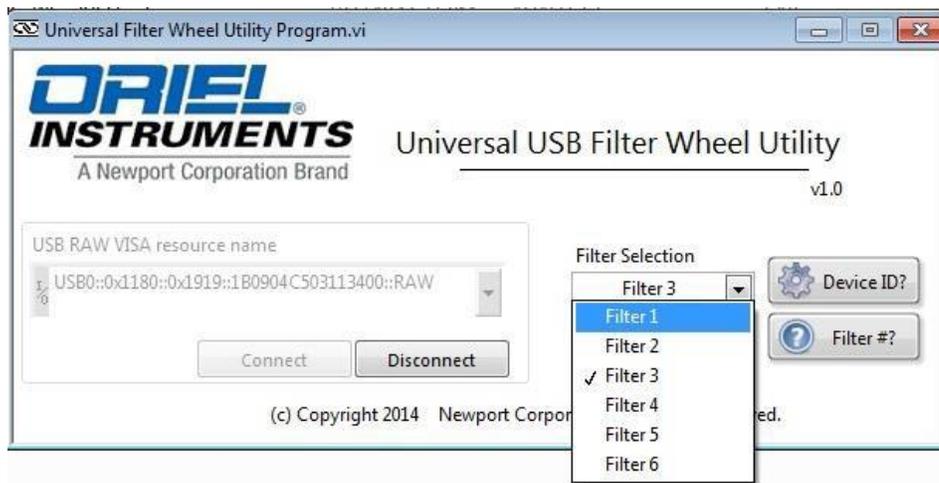


Figure 49: Select Filter

6. To verify which filter is currently in the light path, click “Filter #?”



Figure 50: Query Filter Position

7. To disconnect from communication with the filter wheel, click “Disconnect.”

7.4 COMMAND SET

To ensure latency and guaranteed delivery, all communications with the device uses the control pipeline defined by the USB communications protocol.

COMMAND SYNTAX

Use Line Feed <LF> as the termination character for all commands. The HEX value for <LF> is "0a" or the decimal value is 10

Example:

Command: *IDN?<LF>

Response: APEX2-XE, 001, 1.0<LF>
 where <LF> is the line feed character

COMMAND LIST

All commands listed below automatically assumes the line feed character is used as the termination character.

***IDN?** Identification query. Returns identification string 'Oriel Instruments, APEX2 Illumination System, XXXXX, YYYYY, ZZZZZ' where XXXXX, YYYYY and ZZZZZ are the Model, Serial, and Firmware numbers of the system.

Example:

Command: *IDN?<LF>

Response: APEX2-XE, 001, 1.0

The response indicates that this is a product of Oriel Instruments, the product name is APEX2 Illumination System, the model number is APEX2-XE, the Serial number is 001, and the firmware version is 1.0.

FILT? Queries active filter position. Returns number of the current filter position or -1 for error.

FILT X Sets active filter position, where X is the desired filter position from 1-6.

8 ACCESSORIES

The following is a small list of accessories (sold separately), also available from Oriel Instruments, that are compatible with the APEX2. For more information, visit Newport's website (<http://www.newport.com/>) or contact a Newport Technical Sales Engineer.

8.1 MONOCHROMATOR MOUNTING KITS

Oriel offers a selection of mounting kits which provide a common baseplate and coupling components necessary to connect the APEX2 to many of Oriel's monochromators and spectrographs.

Instrument	Mounting Kit Model
Cornerstone 130(B) Monochromator	APEX2-CS1-KIT
Cornerstone 260 Monochromator	APEX2-CS2-KIT
MS260i Spectrograph	APEX2-CS2-KIT
MS257 Monochromator	APEX2-MS1-KIT
MS257 Spectrograph	APEX2-MS1-KIT

8.2 FIBER FOCUSING ASSEMBLIES

Fiber Focusing Assemblies efficiently collect and focus the collimated beam of light from the APEX2 onto a fiber optic cable or liquid light guide. These assemblies attach directly to the output of the APEX2.



Figure 51: Fiber Focusing Assembly

8.3 FOCUSING LENS ASSEMBLIES

To focus the collimated output of the APEX2 into a concentrated spot, connect the Oriel model 6195 lens holder (shown in Figure 52) to the APEX2 output flange. Choose any 1.5-inch (38.1 mm) lens from Newport's extensive selection or order one of Oriel's preconfigured Focusing Lens Assemblies.



Figure 52: Model 6195 Optic Holder

8.4 BEAM TURNERS

The light path extends horizontally out from the APEX2. To direct the beam up, down or sideways, order one of Oriel's beam turning reflector holders. Then choose a dichroic or full reflector to install into the holder.



Figure 53: Beam Turning Reflector Holder

8.5 UV SAFETY ACCESSORIES

Ultraviolet (UV) light is known to cause both short-term and long-term damage to eyes and skin. The Xenon and QTH lamps utilized in the APEX2 illumination systems emit UV light. Although QTH lamps emit UV light to a lesser degree than Xenon lamps, necessary precautions to protect the eyes and skin from UV exposure should be followed when both QTH and Xenon lamps are being used. Oriel offers various accessories to protect the eyes, face and hands. Figure 54 shows protective gloves, which are available in two different sizes. Figure 55 shows the model 49125 UV Safety Spectacles and 49126 UV Safety Goggles, each capable of providing sufficient UV light protection when operating the APEX2.



Figure 54: Protective Gloves



Figure 55: 49125 and 49126 UV Safety Glasses

9 MAINTENANCE

9.1 LAMP REPLACEMENT

The average life of each lamp type is noted below. Note that the lifetime of the lamp can be affected by the manner in which it is used. Frequent ignitions, contamination of the lamp envelope and an excessively hot operating environment all will lead to reduced lamp life.

To avoid system down time, consider purchasing a replacement lamp as the lamp nears the end of its useful life. Refer to the Section 3.5 of this manual for information on lamp installation.

<u>APEX2 Model</u>	<u>Lamp Type</u>	<u>Lamp Model</u>	<u>Average Lifetime</u>
APEX2-QTH	100 W QTH	6338	500 hours
APEX2-XE	100 W Xe	6252	750 hours

9.2 CLEANING

Clean the exterior of the APEX2 when the illuminator is not hot from use. Use a clean, dry cloth. Ensure that the ventilation holes are not blocked with dust. Vacuum the openings, if necessary, from the outside of the unit. When not in use, cover the output port of the APEX2. Always be sure to remove the output port cover before turning on the lamp.

9.3 LAMP CARE AND HANDLING

Do not allow any contaminants or fingerprints to adhere to the lamp envelope or reflector. Always wear powder-free gloves when handling lamps. Handle the lamp by holding the outside of the reflector only. If a lamp becomes contaminated, do not use it before cleaning its envelope with isopropyl alcohol. Dry completely before using.

If contaminants are not removed, it may lead to reduced light output, overheating, damage to the envelope and premature failure. The lamp's reflector is coated and should not be wiped or touched at all. Remove any particulates such as dust using a bulb blower.

9.4 FUSE REPLACEMENT

The fuse compartment may be accessed from the back of the APEX 2 as shown in Figure 56. Only qualified personnel should access the fuses.



Figure 56: Fuse Compartment

10 SPECIFICATIONS

10.1 APEX2 ILLUMINATOR

Parameter	APEX2-XE Specification	APEX2-QTH Specification
Light Source	100 watt Xenon lamp with integrated reflector	100 watt Quartz Tungsten Halogen lamp with integrated reflector
Lamp Model	6252	6338
Lamp Lifetime (Average)	750 hours	500 hours
External Lamp Intensity Control	None	0 to 5 VDC control signal 0 VDC = no lamp output 5 VDC = 100 watt lamp output
Divergence (Half Angle)	5.0° ± 0.5°	5.0° ± 0.5°
Spot Size (measured at 30 mm from output flange)	41.5 ± 2 mm diameter	41.5 ± 2 mm diameter
Light Ripple (rms)	< 0.2%	< 0.07%
Temporal Stability	< 1%	
Lens Material	Fused Silica	
Lens Transmittance Range	200 to 2500 nm	
Number of Filters	Holds up to six (6) filters, ordered separately. Uses maximum five (5) with MS257 monochromator	
Filter Diameter	1 inch [25.4 mm]	
Filter Thickness	0.23 inch [6.0 mm] maximum	
Connection Flange	Oriell 1.5-Inch Series Female Flange	
Shutter	Manually Controlled	
AC Input and 5x20 mm Fuses (quantity 2)	100 to 120 VAC, 50 to 60 Hz / 1.5 A, 250 V Time Delay Fuse 200 to 240 VAC, 50 to 60 Hz / 0.8 A, 250 V Time Delay Fuse	
Operating Environment	5 to 40°C [41 to 104°F] , indoor use only < 80% relative humidity (non-condensing) < 3000 m [9843 ft] altitude	
IEC Classification	Installation Category II, Pollution Degree 2	
Storage Temperature	0 to 50°C [32 to 122°F]	
Size	22.49 x 6.57 x 8.92 inch [571 x 167 x 227 mm]	
Weight	13 lbs. [5.9 kg]	
Optical Axis Height	4 inch [102 mm]	
Safety Interlock	Lamp shuts off when cover is removed or when internal temperature exceeds normal operating conditions	
CE Compliance	Yes	
RoHS Compliance	Yes	


10.2 UNIVERSAL FILTER WHEEL UTILITY SOFTWARE

Parameter	Specification
Operating System	Microsoft Windows XP, Service Pack 3 Microsoft Windows 7, 32-bit and 64-bit Requires Microsoft .net Framework 4.0
Processor Speed	2 GHz minimum
Random Access Memory (RAM)	1 GB minimum
Peripheral Hardware	USB 2.0 port
Hard Drive Space	800 MB minimum

Note:

One USB 2.0 port is required to connect between the computer and APEX2 filter wheel. Additional USB ports may be required when controlling other instruments, such as a monochromator, lock-in amplifier, power meter, etc.

11 EU DECLARATION OF CONFORMITY

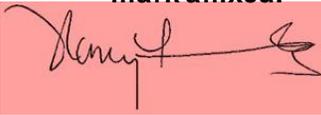EC DECLARATION OF CONFORMITY	
Manufacturer's name:	Newport Corporation
Manufacturer's address:	150 Long Beach Boulevard Stratford, CT 06615 USA
Declares that the product:	
Product Name:	APEX2 Illumination System
Model Numbers:	APEX2-XE, APEX2-QTH
Type of equipment:	Electrical equipment for measurement, control and laboratory use in industrial locations
conforms to the following Product Specifications:	
Safety:	EN 61010-1:2010
EMC:	EN 61326-1:2013
complies with the following Directives:	
	2014/30/EU EMC Directive 2006/95/EC Low Voltage Directive
and accordingly, carries the  mark	
mark affixed:	Beaune; April 10, 2014
	
John Park	Bruno Rety Authorized to compile technical documentation
Engineering Manager, Oriel Products	Group Director, PPT Instrument and Motion Europe
150 Long Beach Boulevard	Micro-Controle Division of Newport Corporation
Stratford, CT 06615 USA	Zone Industrielle

45340 Beaune la Rolande, France

12 WARRANTY AND SERVICE

12.1 CONTACTING ORIEL INSTRUMENTS

Oriel Instruments belongs to Newport Corporation's family of brands. Thanks to a steadfast commitment to quality, innovation, hard work and customer care, Newport is trusted the world over as the complete source for all photonics and laser technology and equipment.

Founded in 1969, Newport is a pioneering single-source solutions provider of laser and photonics components to the leaders in scientific research, life and health sciences, photovoltaics, microelectronics, industrial manufacturing and homeland security markets.

Newport Corporation proudly serves customers across Canada, Europe, Asia and the United States through 9 international subsidiaries and 24 sales offices worldwide. Every year, the Newport Resource catalog is hailed as the premier sourcebook for those in need of advanced technology products and services. It is available by mail request or through Newport's website. The website is where one will find product updates, interactive demonstrations, specification charts and more.

To obtain information regarding sales, technical support or factory service, United States and Canadian customers should contact Oriel Instruments directly.

Oriel Instruments
1791 Deere Avenue
Irvine, CA 92606

Telephone: (800)-222-6440 (toll-free in United States)

Fax: (949) 253-1680

Sales/Technical Assistance: oriel.sales@newport.com

Repair Service: service@newport.com

Customers outside of the United States must contact their regional representative for all sales, technical support and service inquiries. A list of worldwide representatives can be found on Newport's website: <http://www.newport.com/locations/worldwidecontacts.aspx>.

12.2 REQUEST FOR ASSISTANCE / SERVICE

Please have the following information available when requesting assistance or service:

Contact information for the owner of the product.
Instrument model number (located on the product label).
Product serial number and date of manufacture (located on the product label).
Description of the problem.

To help Oriel's Technical Support Representatives diagnose the problem, please note the following:

Is the system used for manufacturing or research and development?
What was the state of the system right before the problem?
Had this problem occurred before? If so, when and how frequently?
Can the system continue to operate with this problem, or is it non-operational?
Were there any differences in the application or environment before the problem occurred?

12.3 REPAIR SERVICE

This section contains information regarding factory service for this product. The user should not attempt any maintenance or service of the system beyond the procedures outlined in this manual. This product contains no user serviceable parts other than what is noted in this manual. Any problem that cannot be resolved should be referred to Oriel Instruments.

If the instrument needs to be returned for service, a Return Material Authorization (RMA) number must be obtained prior to shipment to Oriel Instruments. This RMA number must appear on both the shipping container and the packaged documents.

Return the product to Oriel Instruments, freight prepaid, clearly marked with the RMA number and it either will be repaired or replaced at Oriel's discretion.

Oriel is not responsible for damage occurring in transit. The Owner of the product bears all risk of loss or damage to the returned Products until delivery at Oriel's facility. Oriel is not responsible for product damage once it has left the facility after repair or replacement has been completed.

Oriel is not obligated to accept products returned without an RMA number. Any return shipment received by Oriel without an RMA number may be reshipped by Newport, freight collect, to the Owner of the product.

12.4 NON-WARRANTY REPAIR

For Products returned for repair that are not covered under warranty, Newport's standard repair charges shall be applicable in addition to all shipping expenses. Unless otherwise stated in Newport's repair quote, any such out-of-warranty repairs are warranted for ninety (90) days from date of shipment of the repaired Product.

Oriel will charge an evaluation fee to examine the product and determine the most appropriate course of action. Payment information must be obtained prior to having an RMA number assigned. Customers may use a valid credit card, and those who have an existing account with Newport Corporation may use a purchase order.

When the evaluation had been completed, the owner of the product will be contacted and notified of the final cost to repair or replace the item. If the decision is made to not proceed with the repair, only the evaluation fee will be billed. If authorization to perform the repair or provide a replacement is obtained, the evaluation fee will be applied to the final cost. A revised purchase order must be submitted for the final cost. If paying by credit card, written authorization must be provided that will allow the full repair cost to be charged to the card.

12.5 WARRANTY REPAIR

If there are any defects in material or workmanship or a failure to meet specifications, notify Oriel Instruments promptly, prior to the expiration of the warranty.

Except as otherwise expressly stated in Oriel's quote or in the current operating manual or other written guarantee for any of the Products, Oriel warrants that, for the period of time set forth below with respect to each Product or component type (the "Warranty Period"), the Products sold hereunder will be free from defects in material and workmanship, and will conform to the applicable specifications, under normal use and service when correctly installed and maintained. Oriel shall repair or replace, at Oriel's sole option, any defective or nonconforming Product or part thereof which is returned at Buyer's expense to Oriel facility, provided, that Buyer notifies Oriel in writing promptly after discovery of the defect or nonconformity and within the Warranty Period. Products may only be returned by Buyer when accompanied by a return material authorization number ("RMA number") issued by Oriel, with freight prepaid by Buyer. Oriel shall not be responsible for any damage occurring in transit or obligated to accept Products returned for warranty repair without an RMA number. Buyer bears all risk of loss or damage to the Products until delivery at Oriel's facility. Oriel shall pay for shipment back to Buyer for Products repaired under warranty.

WARRANTY PERIOD

All Products (except consumables such as lamps, filters, etc.) described here are warranted for a period of twelve (12) months from the date of shipment or 3000 hours of operation, whichever comes first.

Lamps, gratings, optical filters and other consumables / spare parts (whether sold as separate Products or constituting components of other Products) are warranted for a period of ninety (90) days from the date of shipment.

WARRANTY EXCLUSIONS

The above warranty does not apply to Products which are (a) repaired, modified or altered by any party other than Oriel; (b) used in conjunction with equipment not provided or authorized by Oriel; (c) subjected to unusual physical, thermal, or electrical stress, improper installation, misuse, abuse, accident or negligence in use, storage, transportation or handling, alteration, or tampering, or (d) considered a consumable item or an item requiring repair or replacement due to normal wear and tear.

DISCLAIMER OF WARRANTIES: EXCLUSIVE REMEDY

THE FOREGOING WARRANTY IS EXCLUSIVE AND IN LIEU OF ALL OTHER WARRANTIES. EXCEPT AS EXPRESSLY PROVIDED HEREIN, ORIEL MAKES NO WARRANTIES, EITHER EXPRESS OR IMPLIED, EITHER IN FACT OR BY OPERATION OF LAW, STATUTORY OR OTHERWISE, REGARDING THE PRODUCTS, SOFTWARE OR SERVICES. NEWPORT EXPRESSLY DISCLAIMS ANY IMPLIED WARRANTIES OF MERCHANTABILITY OR FITNESS FOR A PARTICULAR PURPOSE FOR THE PRODUCTS, SOFTWARE OR SERVICES. THE OBLIGATIONS OF ORIEL SET FORTH IN THIS SECTION SHALL BE ORIEL'S SOLE LIABILITY, AND BUYER'S SOLE REMEDY, FOR BREACH OF THE FOREGOING WARRANTY. Representations and warranties made by any person including distributors, dealers and representatives of Oriel / Newport Corporation which are inconsistent or in conflict with the terms of this warranty shall not be binding on Oriel unless reduced to writing and approved by an expressly an authorized officer of Newport.

12.6 LOANER / DEMO MATERIAL

Persons receiving goods for demonstrations or temporary use or in any manner in which title is not transferred from Newport shall assume full responsibility for any and all damage while in their care, custody and control. If damage occurs, unrelated to the proper and warranted use and performance of the goods, recipient of the goods accepts full responsibility for restoring the goods to their original condition upon delivery, and for assuming all costs and charges.

Confidentiality & Proprietary Rights

Reservation of Title:

The Newport programs and all materials furnished or produced in connection with them ("Related Materials") contain trade secrets of Newport and are for use only in the manner expressly permitted. Newport claims and reserves all rights and benefits afforded under law in the Programs provided by Newport Corporation.

Newport shall retain full ownership of Intellectual Property Rights in and to all development, process, align or assembly technologies developed and other derivative work that may be developed by Newport. Customer shall not challenge, or cause any third party to challenge the rights of Newport.

Preservation of Secrecy and Confidentiality and Restrictions to Access:

Customer shall protect the Newport Programs and Related Materials as trade secrets of Newport, and shall devote its best efforts to ensure that all its personnel protect the Newport Programs as trade secrets of Newport Corporation. Customer shall not at any time disclose Newport's trade secrets to any other person, firm, organization, or employee that does not need (consistent with Customer's right of use hereunder) to obtain access to the Newport Programs and Related Materials. These restrictions shall not apply to information (1) generally known to the public or obtainable from public sources; (2) readily apparent from the keyboard operations visual display, or output reports of the Programs; (3) previously in the possession of Customer or subsequently developed or acquired without reliance on the Newport Programs; or (4) approved by Newport for release without restriction.

First printing 2015

© 2015 by Newport Corporation, Irvine, CA. All rights reserved.
No part of this manual may be reproduced or copied without the prior written approval of Newport Corporation.

This manual has been provided for information only and product specifications are subject to change without notice.
Any change will be reflected in future printings.

Newport Corporation 1791 Deere Avenue Irvine, CA, 92606 USA